# INSTRUCTION MANUAL

## High Current Power Supply

### EPS-200-II

# TABLE OF CONTENTS

	Page
<b>Important User Information</b>	<b>3-4</b>
<b>Section 1    General Information</b>	
1.1    Intended Uses and Specifications	5
1.2    Certification	5
1.3    Set-Up	5-6
1.4    Important Electrical Safety Notice	6
<b>Section 2    Operation</b>	
2.1    Display	7
2.2    Set-Up and Operation	8
<b>Section 3    Maintenance of Equipment</b>	
3.1    Care and Handling	9
3.2    Maintenance	9
<b>Section 4    Troubleshooting &amp; Error Messages</b>	
4.1    Troubleshooting	10
4.2    Error Messages	10-11
<b>Section 5    Equipment</b>	<b>12</b>

## IMPORTANT USER INFORMATION

This Instruction Manual will explain how to use this product safely and effectively. Please read and carefully follow the instruction manual in its entirety.

The triangle/exclamation mark symbol alerts the user of the product to important operational, maintenance, and/or warranty requirements.

The triangle/lightning bolt symbol alerts the user of the product to potentially hazardous electrical exposure.

Failure to adhere to the instructions could result in personal and/or laboratory hazards, as well as invalidate any warranty. Always turn off the DC power source prior to disconnecting power cords from the product. Disconnect power cords from the power source first, and then from the product. For maximum safety, always operate this system in an isolated, low traffic area, not accessible to unauthorized personnel. Never operate damaged or leaking equipment.

## WARRANTY AND LIABILITY

This product was produced utilizing the highest practical standards of materials, workmanship, and design. C.B.S. Scientific warrants that the product has been tested and will meet or exceed published specifications. This warranty is valid only if the product has been operated and maintained according to the instructions provided.

C.B.S. Scientific warrants this product to be free from defects in materials and workmanship under normal service for three years from date of shipment. If the product proves defective during this period, C.B.S. Scientific will repair or replace it at our option, free of charge, if returned to us postage prepaid. This warranty does not cover: damage in transit, damage caused by carelessness, misuse or neglect, normal wear through frequent use, damage caused by solvent corrosion, damage caused by improper handling or user alteration, nor unsatisfactory performance as a result of conditions beyond our control. C.B.S. Scientific shall in no event be liable for incidental nor consequential damages, including without limitation, lost profits, loss of income, loss of business opportunities, loss of use and other related damages, however caused, nor any damage arising from the incorrect use of the product.

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## SECTION 1 General Information

### 1.1 Intended Uses and Specifications

The EPS-200-II Power Supply is intended for use with any electrophoretic device designed to operate below 200 volts and 2000 milliamps DC. Four sets of outputs operate in parallel to provide between 5 and 200 volts in 1 V steps, 0.01 – 2 Amps in 0.01 A steps, and up to 200 watts in the constant voltage mode. Features a 24 hour timer. . A single LED display provides output voltage or current readings.

### 1.2 Certification: EN61010-1-1993 (IEC1010-1), CE

### 1.3 Specifications

Outputs : Voltage --- 5 - 200 VDC adjustable in 1 V steps.

Current --- 10-2000mA DC adjustable in 10mA steps.

Power --- 200 W max.

Input power. Three wire with ground

110V±10% 50/60 Hz or 220V±10% 50/60 Hz

Operation Temperature

0-40°C

Humidity

10% to 90% R.H. Non-condensing

Noise Level

27dB

Regulation:

Ripple:

EPS-200-II

±0.3% @ 200V and 200W

Line Regulation:

EPS-200-II

±0.3 % @ 200V and 200W

From 90 ~ 132VAC or 198 ~ 264 VAC

Load Regulation:

EPS-200-II

±0.5 % 200V for a 50% change in output load

Drift:

EPS-200-II

±0.3 % after 30min warm up at 200V and 200W

Physical Dimensions:

Weight

Net ~2kg (4.4lbs) Gross ~kg (6.6lb)

Length

320mm

Width

240mm

Height

73mm

### 1.4 Set-Up. Power supply is intended for LABORATORY USE ONLY.

(A) Unpacking and Set Up: Unpack and carefully examine the power supply. Report any damage immediately. Do not attempt to operate this device if physical damage is present. Save all packing material if damage is found.

## 1.4 Set-Up---Continued

### (B) Connecting to mains electricity supply

- This apparatus is designed to safety class 1. Before connection to the mains electricity supply, examine the information on the apparatus rating label.
- Ensure that the mains supply is single phase alternating current (a.c.) of the stated frequency (Hz), with Neutral nominally at earth potential.
- Check that the supply voltage is within the stated range. The apparatus rating label states the value of the fuse fitted to the apparatus itself.
- Ensure that the plug or outlet circuit is fitted with an appropriate fuse of higher value.

**WARNING: THIS EQUIPMENT MUST BE EARTHED.**

The wires in the electrical lead are colored in accordance with the following code:

Green	Earth	(Ground)
White	Neutral	(Common)
Black	Live	(Phase)

Connect the wires to a non-reversible plug as follows:

Green wire to the terminal marked E (Earth), G (Ground), or colored Green.

White wire to the terminal marked N (Neutral) or Common.

Black wire to the terminal marked L (Live) or Phase.

### (C) IMPORTANT ELECTRICAL SAFETY NOTES:

Consult a qualified electrician if any doubt or if your supply system has any of the following:

- No Earth
- A Color Code Different from the Above
- Reversible Plugs
- Supply and Return Leads that are Both Above Earth Potential

### (D) Locate:


Place the power supply and cell on a level work surface in an authorized work area. Place the power supply in proximity to the electrophoretic cell(s) with which it is to be connected. Be sure to place the power supply in a safe, dry location with all the controls accessible.

## SECTION 2 Operation

### 2.1 Displays

The front and rear panel keys discussed in Fig-3.1 are used to set up and run the EPS-200-II Power Supply. A beep sound will alarm to inform user when front panel key been pressed. A long beep sound will alarm after the run is complete or if an error occurs. This beep sound will continue to alarm every ten seconds until any key is pressed or the power is turned off.

	<b>KEY</b>	<b>Description</b>
1	ON/OFF	ON/OFF Power Supply
2	CONST	Select voltage or current to be constant
3	MODE	Select the volts, milli amps, or time to be displayed
4	△ ▽	Changes the displayed value of the selected parameter
5	RUN	Starts and Stops the output of power from the Power Supply

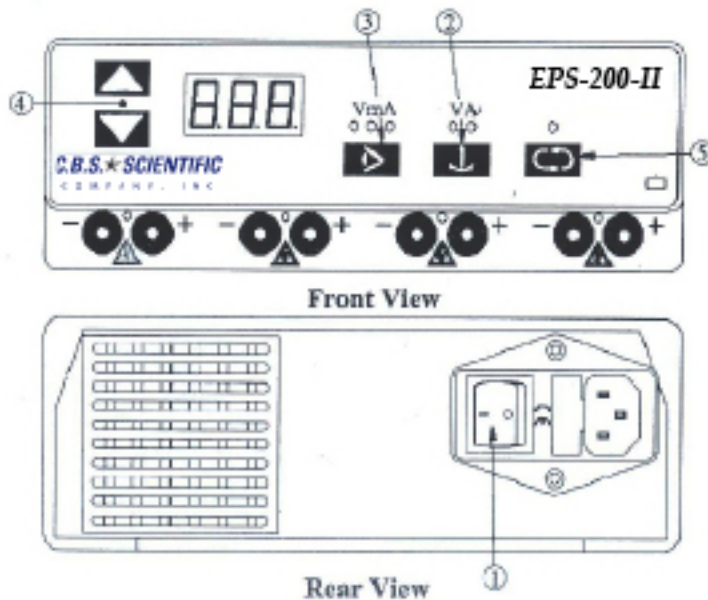1. **MAINS POWER SWITCH**
2. **CONSTANT KEY SELECT VOLTS OR MILLIAMPS**
3. **MODE KEY SELECT.** Display of Volts, Milliamps or Time
4. Changes the displayed value of selected parameter
5. **RUN KEY.** Starts and stops the DC output of power
6. **MAINS POWER CORD RECEPTACLE**

## 2.2 SET UP and OPERATIONS

Step	Action	Description
1	Connect the electrophoresis cell(s) to the power supply.	The power leads are color coordinated to the output terminals.
2	Turn on MAINS power (#1 in Fig. 3.1)	The default display is constant voltage and "0"
3	Select the constant parameter.	V and A are alternately selected by pressing the CONST key (#2 in Fig 3.1).
4	Adjust Limits	Use up or down keys to enter suitable value.
5	Set the default limit if different than maximum default value.	Default current limit is 2000mA Default voltage limit is 200V.
6	If programming a timed run, press MODE to display Time and enter the minutes to be run (countdown timer).	Time is in minutes & the maximum is 24 hours or 1,440 minutes.

CONTINUOUS RUN: Leave the timer at "0" to start a continuous run. The power supply will continue to operate indefinitely, until stopped manually. The timer will start counting up, indicating the length of the run in minutes. After reaching the maximum time of 1440 minutes the timer will automatically reset and recount again from 0. Therefore for over night operation the resulting operation time is the addition of the current timer value with the previous maximum timer 1440 minutes.

TIMED RUN: Press the up and down buttons to set the timer to the desired value (minutes). The timer counts down during the run, indicating the time remaining. At the end of the timed run, a ~3 second beep sounds and the display changes to "OFF". To reset the power supply press any button except the RUN button.

7A POWER FAILURE DETECTION (PFD) The Power Failure Detection feature will save the time, voltage and current settings in the event of a power outage and automatically continue the run when the power is restored. To alert the user that a power failure occurred during the run, the error message "E6" flashes at the end of the run instead of the usual "OFF" message. The "E6" error message means that despite the power outage, the power supply successfully supplied the programmed voltage and current for the entire length of time that was programmed. The time shown at the end of the run is the time the gel ran since power was restored, not the total time of the run. The default settings for Power Failure Detection are:

<u>RUN</u>	<u>Time Settings</u>	<u>PFD</u>	<u>Code</u>
Timed Run	1-1440 minutes	ON	PF1
Continuous Run	0	OFF	PF0

The Power Failure Detection feature can be deactivated during a timed run.

- Press the MODE button to move the indicator light under clock symbol.
- Press the UP and DOWN buttons to enter in the run time in minutes.
- Press both UP and DOWN buttons at the same time. The display will briefly show "PF0", indicating that the PFD mode is disabled.
- Press both the UP and DOWN buttons at the same time again. The display will show "PF1", indicating that the PFD mode is enabled.

8 Press RUN

Start the run.

During a run, you can view the run conditions just press the MODE key. You also can change the value of constant parameter or the length of a timed run by press MODE key and then the up/down keys.  

9 Stop of run

When timer reaches its preset limits the START/STOP key again to manually stop a run. Display will show you OFF.

LIVE PARTS SHOULD NEVER BE EXPOSED UNLESS THE APPARATUS HAS BEEN SWITCHED OFF AND ISOLATED FROM THE MAINS ELECTRICITY SUPPLY.

## **SECTION 3**

### **Maintenance of Equipment**

#### **3.1 Care and Handling**

Use the same precautions as with any electrical device. Do not operate without the cover in place or with any possible short circuit. Do not operate in a damp, humid environment where condensing moisture may short out internal electrical components. When moving to or from a cold room, allow at least eight (8) hours for the unit to temperature equilibrate.

Should this product become excessively warm, smoke, or display erratic information during operation, discontinue use immediately and contact a qualified service representative.

#### **3.2 Maintenance**

The following inspection and maintenance procedures will help maintain the safety and reliable performance. Replacement parts can be ordered by calling 1-858-755-4959 or by contacting your local distributor.

Use only mild, non-abrasive detergents. DO NOT use solvents or immerse in water, as this may cause permanent damage.

- Banana plugs and power cords should be inspected regularly. If the banana plugs become loose or do not feel friction tight replace the plugs or power cords.
- Should power cord assemblies (connectors, wire or shrouds) show any signs of wear or damage (e.g. cracks, nicks, abrasions, or melted insulation), replace them immediately.
- The platinum wire is secured to the banana jack by compression between a stainless washer and the jack nut. The nut/washer interface should be tight and free of corrosion.

## SECTION 4 TROUBLE SHOOTING

### 4.1 Troubleshooting Guide

#### Pre-caution

Wait 10 seconds from time AC power is disconnected to allow high voltage discharge within the power supply.

<b>Symptom</b>	<b>Action and possible causes</b>
Fan is not spinning LED & Display do not light up Front panel power on indicator off	<ol style="list-style-type: none"> <li>1. Check source wall power.</li> <li>2. Change a new, same configure fuse and try again.</li> <li>3. Check if power switch malfunction.</li> <li>4. Contact C.B.S. Scientific Company, Inc.</li> </ol>
Fan is spinning LED & Display do not light up Front panel power on indicator off	<ol style="list-style-type: none"> <li>1. Check source wall power.</li> <li>2. Check if line voltage select error</li> <li>3. Contact C.B.S. Scientific Company, Inc.</li> </ol>
Fan is spinning LED & Display do not light up Front panel power on indicator on	<ol style="list-style-type: none"> <li>1. Contact C.B.S. Scientific Company, Inc.</li> </ol>
Fuse F1 Blown	<ol style="list-style-type: none"> <li>1. Change a new same configure fuse and try again.</li> <li>2. Contact C.B.S. Scientific Company, Inc.</li> </ol>

### 4.2 General Error Messages

<b>Error Code</b>	<b>Explanation</b>	<b>Possible Solution</b>
<b>E1</b>	<b>No load detected</b> The electrophoresis cell(s) are not plugged in. During a run, load current dropped more than 4mA.	<ul style="list-style-type: none"> <li>★ Check all connections</li> <li>★ Check appropriate buffer volume.</li> <li>★ Press the RUN key to resume the run or any other key to clear the error code.</li> </ul>
<b>E2</b>	<b>Unacceptable Value(s) Entered</b> <ul style="list-style-type: none"> <li>★ No value entered</li> <li>★ Voltage below 10V</li> <li>★ Current below 4mA</li> </ul>	Enter acceptable values and press run key to resume the run or any other key to clear the error code.
<b>E3</b>	<b>Change in Load Resister</b> <ul style="list-style-type: none"> <li>★ Electrophoresis cell(s) were added or removed during a run.</li> <li>★ Buffer leakage</li> <li>★ Loose connection in a connected cell.</li> </ul>	Check and correct any potential resistance problems Then press RUN key to resume the run or any other key to clear the error code.


## 4.2 General Error Messages--Continued

<u>Error Code</u>	<u>Explanation</u>	<u>Possible Solution</u>
E4	<b>Power Failure During a Non-Timed Run</b>	Press any key (other than RUN) to clear the error code. Then press the RUN Key to resume the run.
E5	<b>Power Failure During A Timed Run (PFD OFF)</b>	Press any key (other than RUN) to clear the error code. Then press the RUN key to restart the run.
E6	<b>Power Failure During A Timed Run (PFD ON)</b>	Press any key (other than RUN) to clear the error code.
E7	<b>Short Circuit</b> Load current exceeded 500mA	<ul style="list-style-type: none"> <li>★ Check for any short circuit.</li> <li>★ High buffer concentration</li> <li>★ Press the RUN key to resume the run or any other key to clear the error code.</li> </ul>
CHC	<b>Current differential alarm</b> Difference current value (>100mA) between start and final run.	★ Check buffer volume
<b>Any Other Error Code</b>		Contact C.B.S. Scientific Company, Inc.

## 4.3 General Error Messages--Continued

### Replacing the fuses


1. Unplug the power cord from the power supply.
2. Locate the fuse compartment on the rear panel.
3. Use a screwdriver to pry out the fuse compartment. There are two fuses in the compartment. Check and replace any blown fuses with a fuse of the same size and ratings.
4. FUSES (2); 250 V, 5 A, 5 x 20 mm fast blow.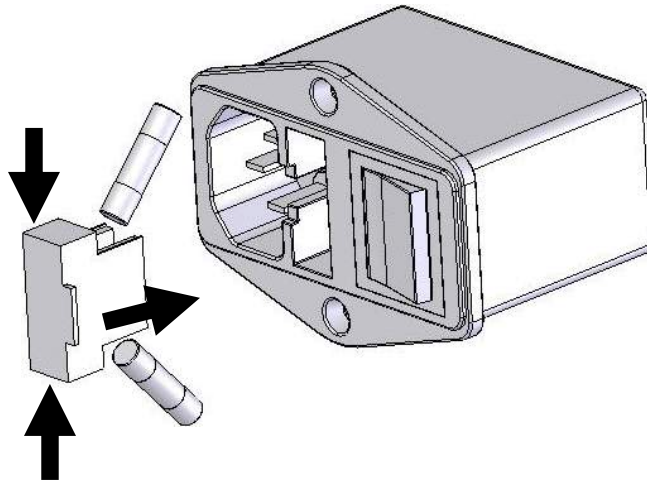

**SECTION 5**  
**Power Supply**

Cat. #	Item
EPS-200-II	Mini Power Supply, 10-300 Volts, 500mA, 110V
EPS-200-IIV	Mini Power Supply, 10-300 Volts, 500mA, 220V

Section 7 — Circuit Diagram  
As special request.

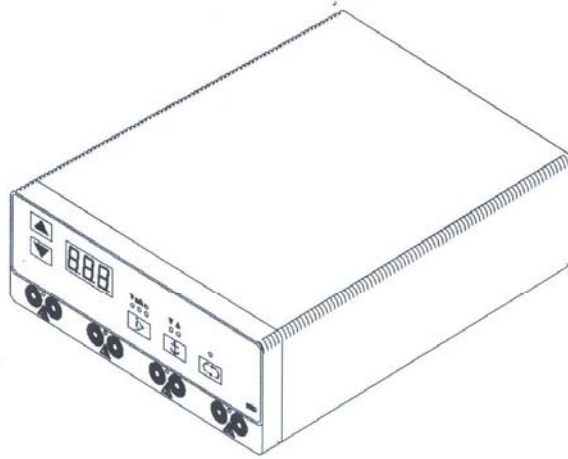# CONTACT INFORMATION

**Telephone:**  
Local or International  
858-755-4959  
Toll Free: 800-243-4959

**Fax:** 858-755-0733

**Online ordering:**  
[www.cbsscientific.com](http://www.cbsscientific.com)

**E-mail address:**  
[sales@cbssci.com](mailto:sales@cbssci.com)

**Mailing address:**  
C.B.S. Scientific Company  
P.O. Box 856  
Del Mar, CA 92014

**Shipping address:**  
C.B.S. Scientific Company  
420 South Cedros  
Solana Beach, CA 92075

**Credit Card Options:**  
Visa/Mastercard  
Discover/American Express