INSTRUCTION MANUAL

Electro-Eluter/Concentrator

ECU-040-20

ECU-040-50

ECU-040-25

ECU-040

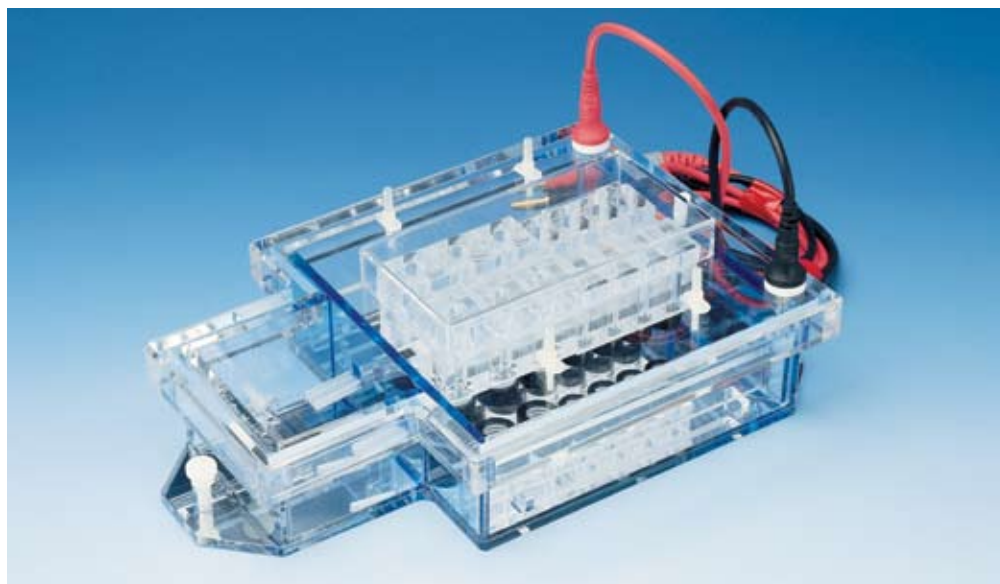
TABLE OF CONTENTS

	Page
Important User Information	3-4
Section 1 General Information	
1.1 Introduction	5
1.2 Specifications	5
1.3 Safety	6
Section 2 Description of parts	
2.1 Unpacking	7
2.2 Components and Assembly	7
Section 3 Instructions for Use	
3.1 Unit Preparation	8
3.2 Elution Block Set-up	8
3.3 Procedure for Protein Elution	8-10
3.4 Procedure for Elution of DNA	11
Section 4 Maintenance of Equipment	
4.1 Care and Handling	12
4.2 Maintenance	12
Section 5 Ordering Information	13

IMPORTANT USER INFORMATION

This Instruction Manual will explain how to use this product safely and effectively. Please read and carefully follow the instruction manual in its entirety.

The triangle/exclamation mark symbol alerts the user of the product to important operational, maintenance, and/or warranty requirements.

The triangle/lightning bolt symbol alerts the user of the product to potentially hazardous electrical exposure.

Failure to adhere to the instructions could result in personal and/or laboratory hazards, as well as invalidate any warranty. Always turn off the DC power source prior to disconnecting power cords from the product. Disconnect power cords from the power source first and then from the product. For maximum safety, always operate this system in an isolated, low traffic area, not accessible to unauthorized personnel. Never operate damaged or leaking equipment.

WARRANTY AND LIABILITY

This product was produced utilizing the highest practical standards of materials, workmanship, and design. C.B.S. Scientific warrants that the product has been tested and will meet or exceed published specifications. This warranty is valid only if the product has been operated and maintained according to the instructions provided.

C.B.S. Scientific warrants this product to be free from defects in materials and workmanship under normal service for one year from date of shipment. If the product proves defective during this period, C.B.S. Scientific will repair or replace it at our option, free of charge, if returned to us postage prepaid. This warranty does not cover: damage in transit, damage caused by carelessness, misuse or neglect, normal wear through frequent use, damage caused by solvent corrosion, damage caused by improper handling or user alteration, nor unsatisfactory performance as a result of conditions beyond our control. C.B.S. Scientific shall in no event be liable for incidental nor consequential damages, including without limitation, lost profits, loss of income, loss of business opportunities, loss of use and other related damages, however caused, nor any damage arising from the incorrect use of the product.

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SECTION 1 General Information

1.1 Introduction

The Electro-Eluter/Concentrator is designed to elute and concentrate macromolecules simultaneously from acrylamide or agarose gels. Typical applications include recovery of proteins from SDS-acrylamide gels, and of restriction enzyme-treated DNA from agarose gels. The elution of proteins from gel strips has become a useful technique for preparing proteins for primary structure analysis because this procedure can be used after Coomassie staining to visualize the location of the protein bands in the gels.

Eluter/Concentrator blocks consist of a precision-machined sample recovery channel and two silicone-Teflon® washer/open-top cap assemblies for holding dialysis membranes securely in place.

The three-chamber Electro-Eluter is designed to allow buffer cycling. The buffer drain outlets on the side for the two electrode chambers allow overflow buffer to drop into the mixing chamber. The mixed buffer is then pumped back into the electrode chambers. This prevents build-up of a pH gradient, which can occur during long-term electrophoresis, interfering with the ionic migration of the protein or DNA sample. The buffer cycling system also eliminates the need to change buffer during electrophoresis.

1.2 Specifications

Constructions

Buffer chamber, lid	Acrylic, mirror
Electrodes	Platinum wire .012" diameter
Power cords	FR Urethane, rated 7500V, 200mA, 65° C
Blocks	Acrylic
Caps	Bakelite*
Washers	TFE/Silicone
Safety Certification	EN61010-1-1993 (IEC1010-1)

	ECU-040	ECU-040-25	ECU-040-20	ECU-040-50	ECB-102	ECB-105
Shipping weight	7lbs	8lbs	8lbs	8lbs	2lbs	2lbs
Overall Size (l)x(w)x(h) cm	30.8x20.3x12.7	30.8x20.3x12.7	30.8x20.3x12.7	30.8x20.3x12.7	1.9x5.7x2.54	5.7x5.7x2.54
Recommended buffer volume:	550	550	550	550	-	-
Loading chamber capacity	—	—	—	—	2.0ml	5.0ml
Migration channel volume	—	—	—	—	5.0ml	12.0ml
Sample recovery volume	—	—	—	—	200ul	500ul
Distance between electrodes	15cm	15cm	15cm	15cm	—	—
Voltage limit	250	250	250	250	—	—

1.3 Safety

Power to the Electro-Eluter/Concentrator unit is to be supplied by an external DC voltage power supply that must be ground isolated so that the DC voltage output floats with respect to ground. For any power supply used, the maximum specified operating parameters for the units are:

Maximum Limits

250 VDC voltage

25 watts power

50 mA current

60° C ambient temperature

Current to the unit, provided from the external power supply, must enter the unit through the lid assembly, providing a safety interlock to the user. Current to the unit is broken when the lid is removed. **Do not attempt to use the unit without the safety lid. Always turn the power supply off before removing the lid, or when working with the unit in any way. Follow safety precautions specified by the power supply manufacturer.**

Description of Parts


2.1 Unpacking

Please verify that your Electro-Eluter Concentrator comes complete with the following components:

- Buffer Chamber with leveling base, spirit level, platinum electrodes, and 2 safety/vapor covers
- Power leads
- 3 “Y” and 4 straight tubing connectors
- 4 Elution/Concentrator Chambers or “Blocks”. Each Block has 2 black open-top screw caps, 2 Teflon-Silicone flat washers.

Notes: We do not supply dialysis membrane. Use non-charged, well washed membrane such as Spectrapore. A small peristaltic pump must be used for buffer re-circulation.

Electro-Eluter Components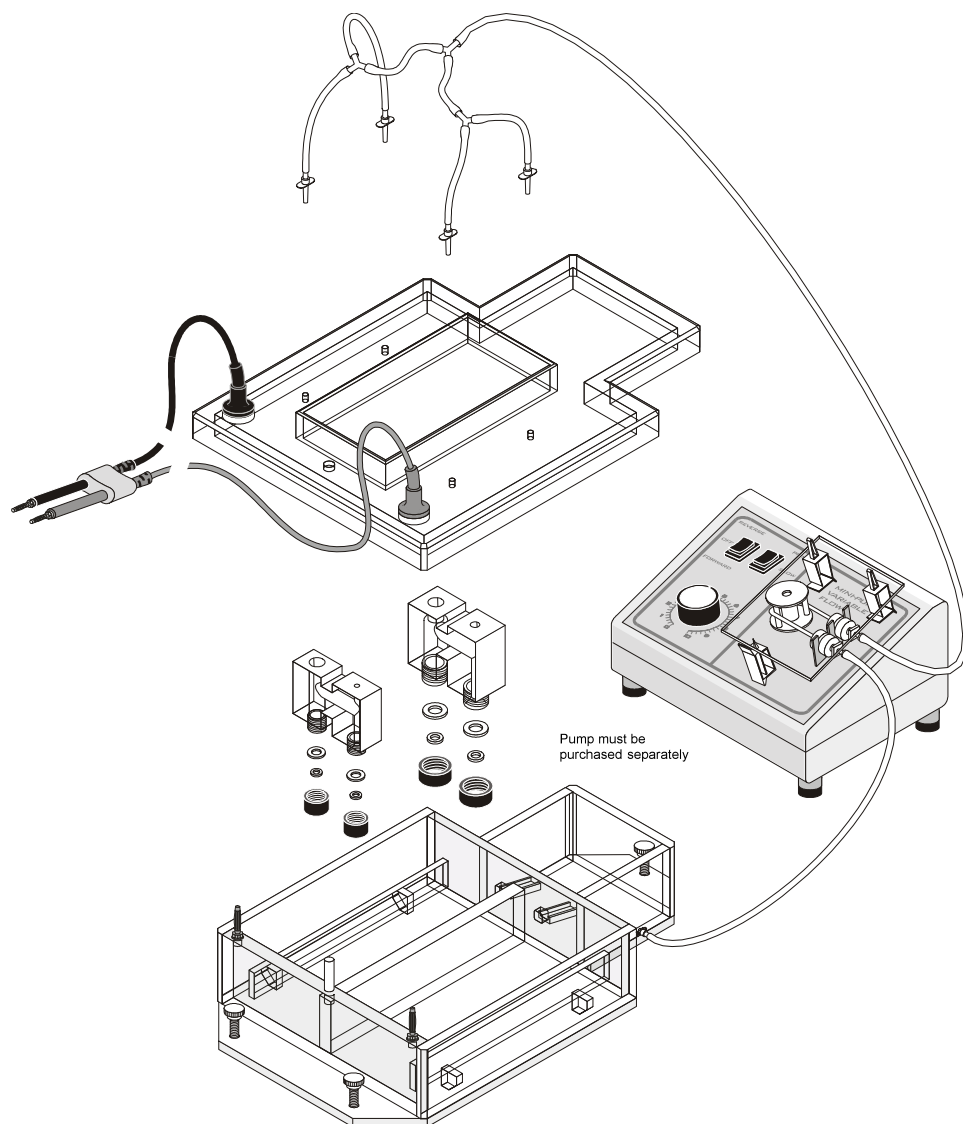

SECTION 3


Instructions for Use

3.1 Unit Preparation

Place the Electro-Eluter Concentrator and the Elution blocks on a level work surface in an authorized work area.

3.2 Elution Block Set-Up

To assemble Block for electroelution, unscrew black caps and remove washers. Cut dialysis membrane into circles the same diameter as the Teflon/Silicone Washers. Place the membrane onto the slightly tan side of the washer inside the black cap (the Teflon will allow the membrane to slide instead of binding on the washer). Screw the black cap with the washer and membrane into place as shown in figure 1.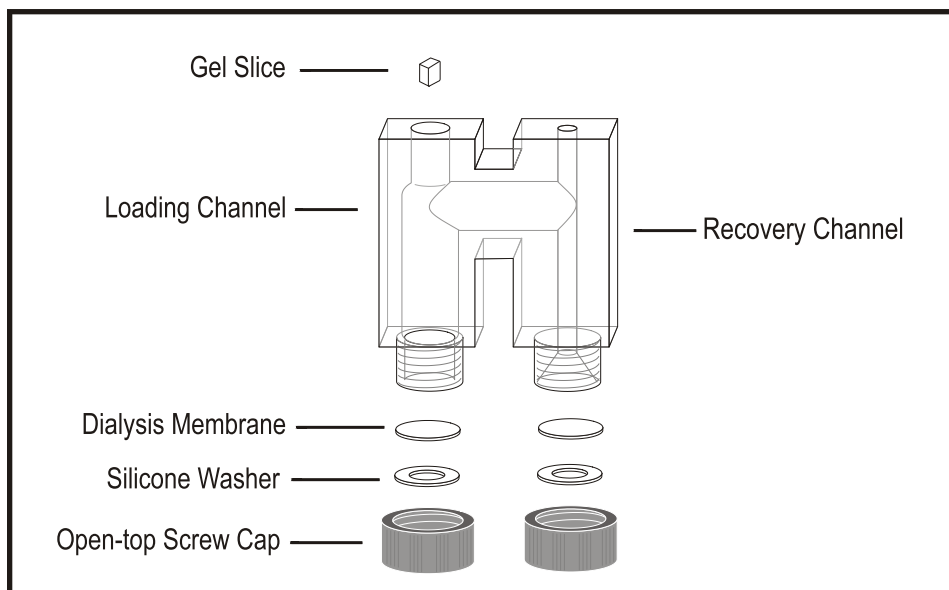


3.3 Procedure for Protein Elution


1. For electroelution of protein samples prior to sequencing, read carefully the Hunkapillar et. al. paper in Methods in Enzymology, Vol. 91, 1983, pgs. 227-236.
2. Equilibrate each block with 1% Ethanolamine to take away most of the non-specific binding sites for protein attachment. The Plexiglas walls of the blocks, if not treated, may decrease the yield from electroelution by non-specific absorption. Dilute the Ethanolamine in running buffer.
3. The gel slice should be equilibrated in running buffer, as per the Hunkapillar paper. On the large volume side of the Block, top load the gel slice onto the membrane. Fill the entire chamber with running buffer $\frac{3}{4}$ of the way up to the cross channel. Tap blocks with finger so that ALL THE AIR BUBBLES TRAPPED INSIDE ARE DISLODGED.
4. Insert loaded blocks into unit. Flush air trapped beneath each black cap with syringe and bent needle. Be careful not to puncture membrane.

5. **Buffer Cycling System Set-up.** This unit is designed to maximize yield by preventing any pH gradient from forming inside the block by constantly re-cycling buffer. Attach tubing to outlet on side of the third chamber to dual peristaltic pump. Split outlets from pump with “Y” connectors, followed by attachment to the tubing connectors mounted in the safety covers. The flow of recycled buffer drops out of anode and cathode buffer chambers into the mixing chamber, and is then returned via the pump system. The remixed buffer drops onto the small plastic strips, which protect the platinum wire. This prevents unwanted small air bubbles from re-entering the buffer.

Buffer Re-Circulation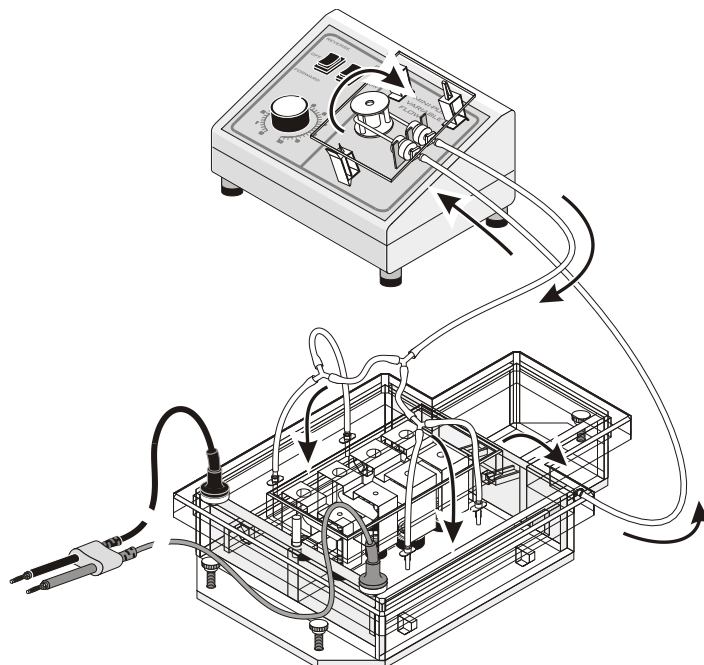

6. Running conditions for
 - a) **Proteins prepared for Sequencing: Electrophoresis** is carried out as per Hunkapillar et. al.
 - b) **Elution of Protein from SDS-Acrylamide Gels, not to be Sequenced:** Buffer: 50mM Tris Acetate (pH 7.8), 1mM EDTA, Current: 60-70 Volts, Duration: 15-20 hours
7. At the end of electrophoresis remove the blocks carefully from the chamber. Pipette off and discard most of the running buffer in the loading and cross channels. Re-suspend the eluted sample on the node membrane, typically 100-200ul of buffer (500ul in the 5ml block). For initial runs save all of the buffer in the collection side above the conical chamber. If the elution is incomplete, protein sample may not have migrated all the way down the membrane. Therefore, check yield before discarding collection side buffer. **See diagram on page 10.**

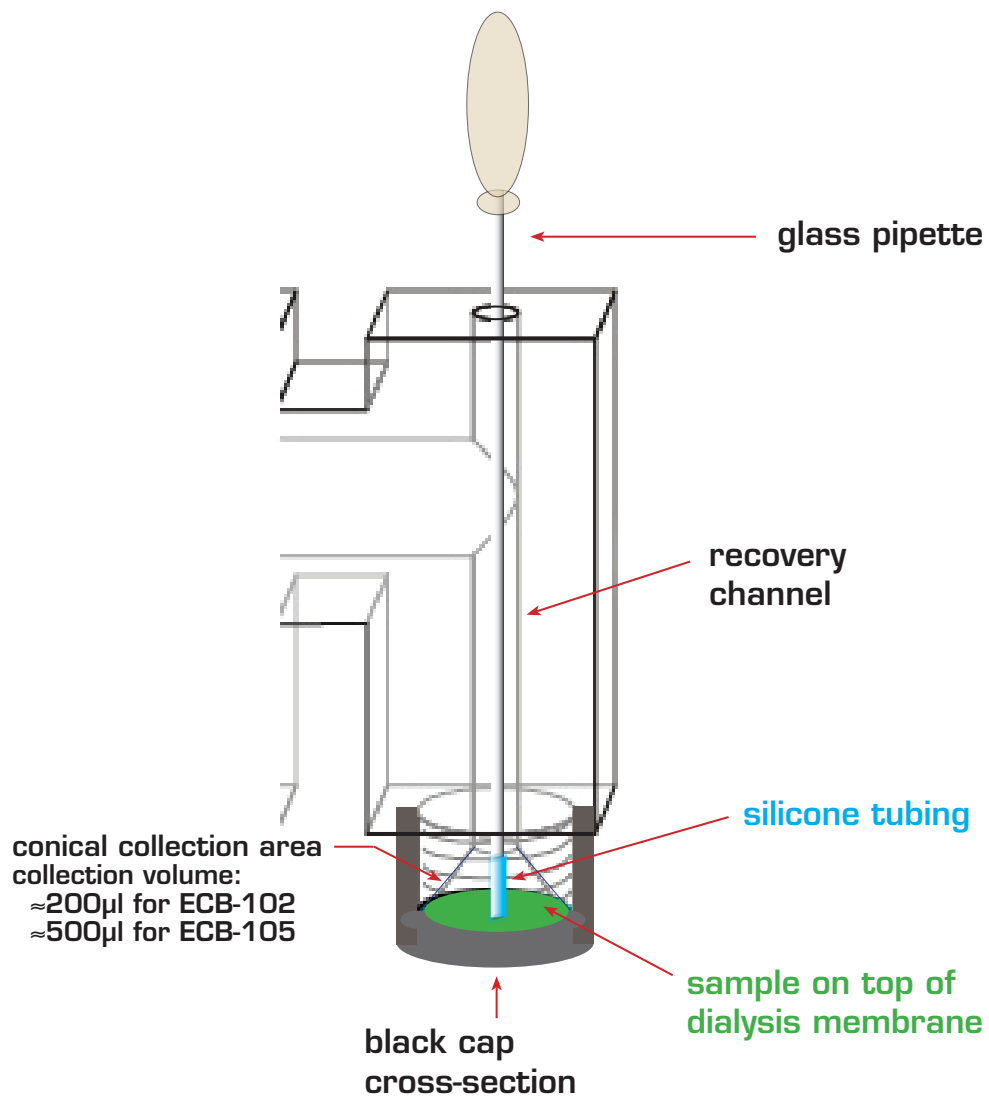3.4 Procedure for Elution of DNA

1. Please read through directions for elution of proteins. Follow instructions 3-5 of Section 3.1 using the buffer suggestions below.
 - a. Buffer: 5mM Tris-HoAc pH 7.5, 1mM EDTA, Current: 100-150 Volts @ 10mA, Duration: Minimum of 5 hours for large DNA (8-10 kbp); 1.5 hours for 300 bp.
 - b. Comparisons of recoveries after 5 hours to 18 hours are similar. Therefore, it does not hurt to run overnight when running large DNA fragments.
2. All electroelution preps are extensively Phenol extracted after run (2x, 1 hour) followed by chloroform ether extraction. This is to remove any contaminants including dissolved agarose. **Note:** Most people use a Brom-phenol-Blue Sucrose marker in with the piece of agarose to make sure current flow is unobstructed by bubbles. After loading sample into chamber, drop a little of this marker into the bottom of the loading chamber. Watch for movement of dye after power is turned on.
3. DNA recovered from our ECU units has been tested for a number of viable characteristics including kinase treatment, re-restriction, and ability to transform E. Coli and has been found to be more active than DNA recovered by conventional techniques. Also, the percent recoveries of DNA and the percent of specific E. Coli transformation by that DNA are much better than traditional methods, despite extensive Phenol extraction.

SECTION 4 Maintenance of Equipment

4.1 Care and Handling

The plastic components of the Electro-Eluter Concentrators are fabricated from acrylic. Electrodes and connectors are made from pure platinum, stainless steel, and chrome plated brass. As with any laboratory instrument, adequate care ensures consistent and reliable performance.

After each use, rinse buffer chamber with de-ionized water. Wipe dry with a soft cloth or paper towel, or allow to air dry. Whenever necessary, all components may be washed gently with water and a non-abrasive detergent, and rinsed and dried as above. *Never* use abrasive cleaners, glass cleaning sprays or scouring pads to clean the components, as these will damage the unit and components.

Additional precautions:

- Do not autoclave or dry-heat sterilize the apparatus or components.
- Do not expose the apparatus or components to phenol, acetone, benzene, halogenated hydrocarbon solvents or alcohols.
- Avoid prolonged exposure of the apparatus or components to UV light.
- Do NOT treat with diethylpyrocarbonate (DEPC)-treated water for extended periods at 37°C. A brief rinse with DEPC-water is sufficient after a thorough wash.

4.2 Maintenance

The following inspection and maintenance procedures will help maintain the safety and reliable performance of the Electro-Eluter Concentrator. Replacement parts can be ordered by calling 1-858-755-4959 or by contacting your local distributor.

- Banana plugs and power cords should be inspected regularly. If the banana plugs become loose or do not feel friction tight replace the plugs or power cords.
- Should power cord assemblies (connectors, wire or shrouds) show any signs of wear or damage (e.g. cracks, nicks, abrasions, or melted insulation), replace them immediately.
- The platinum wire is secured to the banana jack by compression between a stainless washer and the jack nut. The nut/washer interface should be tight and free of corrosion.

SECTION 6 Equipment and Accessories

<u>Cat. #</u>	<u>Item</u>
ECU-040-20	Electro-Eluter Concentrator. Includes four 2.0ml blocks
ECU-040-50	Electro-Eluter Concentrator. Includes four 5.0ml blocks
ECU-040-25	Electro-Eluter Concentrator. Includes two 2.0ml blocks and two 5.0ml blocks
ECU-040	Electro-Eluter Concentrator. Unit only
ECB-102	Additional 2.0ml Eluter/Concentrator block
ECB-105	Additional 5.0ml Eluter/Concentrator block
ECB-102-C	Additional 2.0ml Cap
ECB-105-C	Additional 5.0ml Cap
ECB-102-W	Additional 2.0ml Washer
ECB-105-W	Additional 5.0ml Washer

Power Supply

<u>Cat. #</u>	<u>Item</u>
EPS-200-II	High Current Mini Power Supply with timer, CE, CV or CC 5-200 Volts, 110V/50-60Hz, current range: 2000mA, 200 Watts
EPS-200-IV	High Current Mini Power Supply with timer, CE, CV or CC 5-200 Volts, 220V/50-60Hz, current range: 2000mA, 200 Watts
EPS-300-II	Mini Power Supply with timer, CE, CV or CC 10-300 Volts, 110V/50-60Hz, current range: 4-500mA, 90 Watts
EPS-300-IV	Mini Power Supply with timer, CE, CV or CC 10-300 Volts, 220V/50-60Hz, current range: 4-500mA, 90 Watts

Casting Pump

<u>Cat. #</u>	<u>Item</u>
MPP-100	Mini Casting/Recirculation Pump

NOTES

NOTES

CONTACT US

Telephone:
Local or International
858-755-4959
Toll Free: 858-755-0733

Fax: 858-755-0733

Check out our new and improved online ordering!

www.cbsscientific.com

E-mail address:
sales@cbssci.com

Mailing address:
C.B.S. Scientific Company
P.O. Box 856
Del Mar, CA 92014

Shipping address:
C.B.S. Scientific Company
10805 Vista Sorrento Pkwy
Suite 100
San Diego, CA 92121

Credit Card Options:
Visa/Mastercard
Discover/American Express