INSTRUCTION MANUAL

Horizontal Midi-Gel Productivity Package

GCMGU-402T

GCMGU-502T

GCMGU-602T

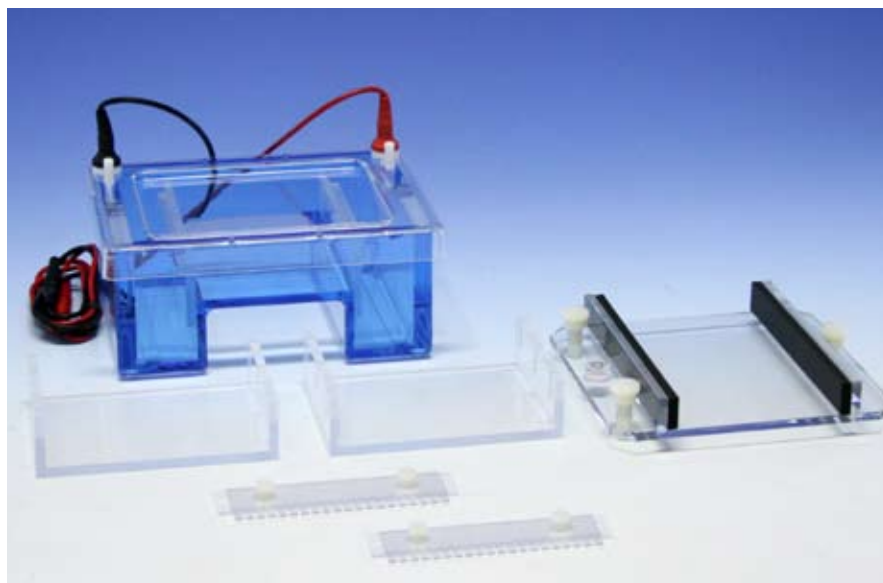
TABLE OF CONTENTS

| | Page |
|--|-------|
| Important User Information | 3-4 |
| Section 1 General Information | |
| 1.1 Introduction | 5 |
| 1.2 Specifications | 5 |
| 1.3 Safety | 6 |
| Section 2 Description of parts | |
| 2.1 Unpacking | 6 |
| 2.2 Components/Assembly | 6 |
| Section 3 Instructions for Use | |
| 3.1 Midi-Gel Unit Preparation and Gel Casting | 7-8 |
| 3.2 Running the Gel | 8 |
| 3.3 Removing the Gel | 9 |
| 3.4 Maximum Well/Comb volumes | 10 |
| Section 4 Running Conditions | |
| 4.1 Recommended Power | 10 |
| 4.2 Recommended Buffers | 10-11 |
| Section 5 Maintenance of Equipment | |
| 5.1 Care and Handling | 11 |
| 5.2 Maintenance | 11 |
| Section 6 Equipment and Accessories | 12-13 |

IMPORTANT USER INFORMATION

This Instruction Manual will explain how to use this product safely and effectively. Please read and carefully follow the instruction manual in its entirety.

The triangle/exclamation mark symbol alerts the user of the product to important operational, maintenance, and/or warranty requirements.

The triangle/lightning bolt symbol alerts the user of the product to potentially hazardous electrical exposure.

Failure to adhere to the instructions could result in personal and/or laboratory hazards, as well as invalidate any warranty. Always turn off the DC power source prior to disconnecting power cords from the product. Disconnect power cords from the power source first, and then from the product. For maximum safety, always operate this system in an isolated, low traffic area, not accessible to unauthorized personnel. Never operate damaged or leaking equipment.

WARRANTY AND LIABILITY

This product was produced utilizing the highest practical standards of materials, workmanship, and design. C.B.S. Scientific warrants that the product has been tested and will meet or exceed published specifications. This warranty is valid only if the product has been operated and maintained according to the instructions provided.

C.B.S. Scientific warrants this product to be free from defects in materials and workmanship under normal service for one year from date of shipment. If the product proves defective during this period, C.B.S. Scientific will repair or replace it at our option, free of charge, if returned to us postage prepaid. This warranty does not cover: damage in transit, damage caused by carelessness, misuse or neglect, normal wear through frequent use, damage caused by solvent corrosion, damage caused by improper handling or user alteration, nor unsatisfactory performance as a result of conditions beyond our control. C.B.S. Scientific shall in no event be liable for incidental nor consequential damages, including without limitation, lost profits, loss of income, loss of business opportunities, loss of use and other related damages, however caused, nor any damage arising from the incorrect use of the product.

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General Information

1.1 Introduction

The mid-range Horizontal Midi-Gel units, MGU-402T (11 x 11cm), MGU-502T (10 x 14cm) and MGU-602T (14 x 14cm) effectively allow rapid separation and identification of various fragments from DNA, RNA, PCR™ products and synthetic oligonucleotides. When used in combination with two or more sample combs, designed for multi-channel pipette loading, multiple screening of large sample numbers can be facilitated easily. The unique tapered baffle system allows immediate gel casting directly onto removable trays without having to seal with agarose plugs or tape. The extensive variety of comb formats and material types ensure a fast and easy method of identifying critically important samples while using minimal quantities of expensive agarose.

Two methods for gel formation: Option 1, in Situ tapered baffle systems for simplified gel casting. The tapered baffle system allows gel casting directly onto the unit or removable trays without having to seal with agarose plugs or tape. Option 2: External horizontal gel casters. Precast your agarose gels for future use in an easy to use and easy to store fashion. These support trays have leveling bases and are lined with soft, coated neoprene to provide firm yet supple support for casting gels without leakage.

1.2 Specifications

Constructions:

| | |
|--------------------------------|---|
| Buffer chamber | Acrylic |
| Safety Cover | Acrylic |
| Electrodes | Platinum wire .012" diameter |
| Power cords | FR Urethane rated 7500VDC, 200mA, 65°C |
| Tapered baffles | Acrylic |
| Combs | Polycarbonate or Teflon |
| Adjustable comb backing | Acrylic |
| UVT Gel Tray | UV transparent acrylic |
| Safety Certification | EN61010-1-1993 (IEC1010-1) |

| | MGU-402T | MGU-502T | MGU-602T |
|---|----------------|------------------|------------------|
| Shipping weight | 4lb. | 6lb. | 7lb. |
| Overall Size (l)x(w)x(h) cm | 22.5 x 13 x 14 | 21.5 x 16.7 x 14 | 27.5 x 16.7 x 14 |
| Gel Size cm | 11 x 11 | 14 x 10 | 14 x 14 |
| Recommended buffer volume: with gel tray | 550 ml | 700 ml | 750 ml |
| Distance between electrodes | 20 cm | 19 cm | 25 cm |
| Voltage limit | 250VDC | 250VDC | 250VDC |

1.3 Safety

Power to the MGU-402T, MGU-502T or MGU-602T unit is to be supplied by an external DC voltage power supply that must be ground isolated, so that the DC voltage output floats with respect to ground. For any power supply used, the maximum specified operating parameters for the units are:

Maximum Limits

250VDC voltage

15 watts power

60 mA current

55°C ambient temperature

Current to the unit, provided from the external power supply, must enter the unit through the lid assembly, providing a safety interlock to the user. Current to the unit is broken when the lid is removed. **Do not attempt to use the unit without the safety lid, and always turn the power supply off before removing the lid, or when working with the unit in any way. Follow safety precautions specified by the power supply manufacturer.**

SECTION 2


Description of Parts

2.1 Unpacking

Please verify that your unit comes complete with the following components:

- Midi-Horizontal electrophoresis unit
- Tapered baffles (2)
- Safety cover with attached DC power leads
- UV Transparent slotted gel tray (2)
- Two combs with adjustable height backing
- One gel casting tray, single place (not pictured)

2.2 Components/Assembly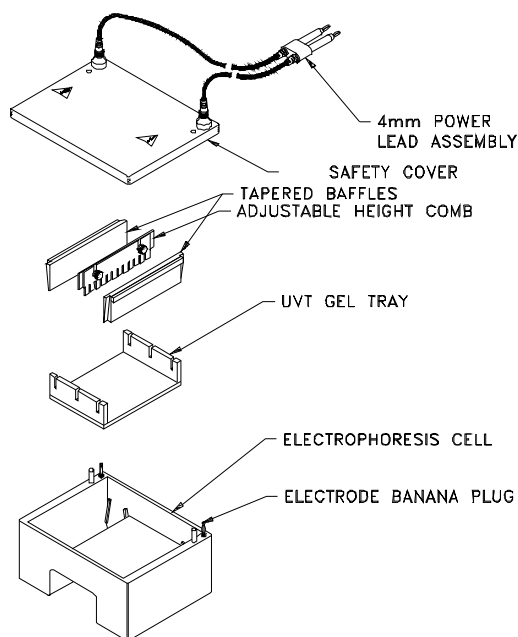

SECTION 3 Instructions for Use

3.1 Midi-Gel Unit Preparation and Gel Casting



Place Midi-Gel unit on a level work surface in an authorized work area. Place the adjustable height comb assembly in the gel tray. Loosen the white screws holding the comb to the backing and adjust comb depth (generally 2mm). Gently tighten the adjusting screws and set the comb aside. **See Table 1 for comb volumes.**


The Midi-Gel unit provides the user with two options for gel casting: (a) **gel casting directly into unit using gel tray** or (b) **gel casting in gel tray outside of unit.**

A. Casting Directly into Unit Using Gel Tray.


1. Place gel tray onto raised bed of unit. Insert the tapered baffles, simultaneously on either side of gel tray and slide into the grooves next to the bed, handles to the outside. The tapered baffles will rest friction tight against the gel tray. **DO NOT** press baffles too tightly into stops.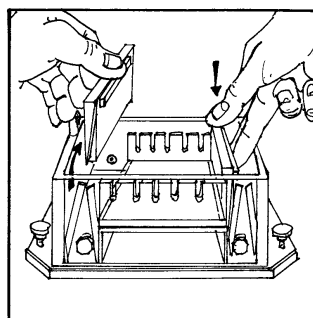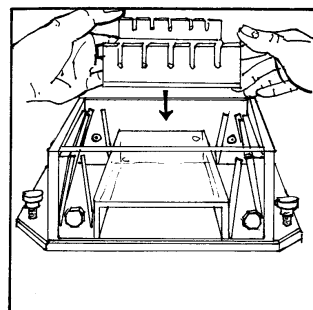

2. Position the gel comb(s). Remembering, current travels from negative (black cathode) towards positive (red anode). Pour the agarose solution onto the gel bed. Allow the agarose slab to cool uncovered until it turns opaque. Agarose around combs will take a few more minutes to cool than other parts of the slab.

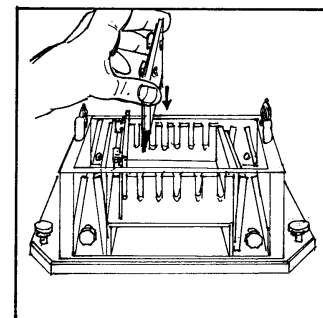
3. Slide the baffles out of their slots and store in the slots provided at either end of the unit near both terminals.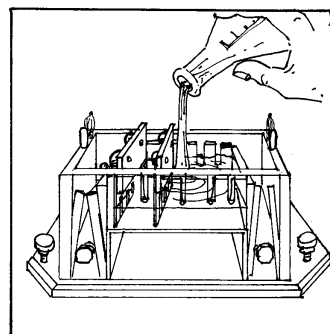


B. Gel Casting into Gel Tray Outside of Unit.

1. Tape both ends of gel tray with tape.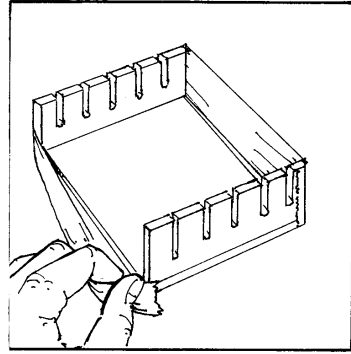

2. Place gel comb(s) into desired slot(s) in the gel tray. Using a pipette or a pipettor, run a bead of agarose solution along the edges of the gel tray where it meets the tape to assure a leak-proof seal. Then pour the agarose solution onto the gel tray. Allow the agarose slab to cool uncovered until it turns opaque. NOTE: (Optional storage) Pre-cast gels can be stored for up to one month if wrapped in plastic wrap and stored at 4°C. Remove combs prior to storing.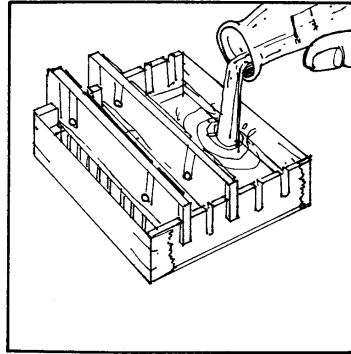

3. Remove gel tape at each end. Position the tray in the unit, remembering, current travels from negative (black cathode) towards positive (red anode).


C. Gel Casting with External Gel Caster.

1. Place external gel caster on level surface and level using thumb screws.
2. Push gel tray against one side of gel caster, then the other. Confirm it the gel tray is flat to the bottoms of the caster


3.2 Running the gel


1. Add enough buffer to fill both reservoirs and overflow the surface of the gel to a depth of 2-3mm. Gently remove comb(s). Flush out any air bubbles in the wells. Load the samples into the sample wells. NOTE: DO NOT FORGET TO LOAD DNA SIZE STANDARD.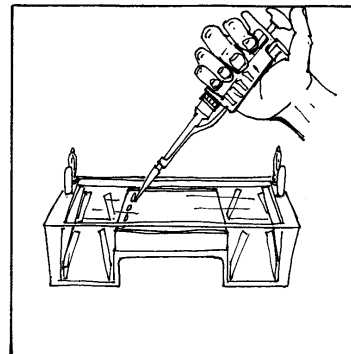

- Align safety cover over the unit and carefully attach, so not to disturb samples.

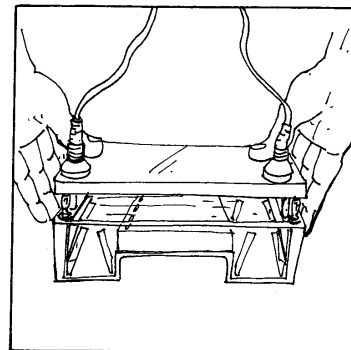
- Connect the leads to the power supply, matching the color-coded red to red and black to black. **See Section 4.1 for recommended power conditions.** Begin separation by electrophoresis.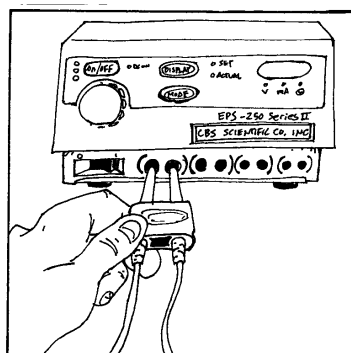

3.3 Removing the Gel

- Turn the power supply off and disconnect the leads from the power supply.
- Remove the safety cover from the unit, by placing thumbs on white posts next to red & black connectors, then pushing down while pulling up with fingers under lid. **DO NOT pull on power cords.**

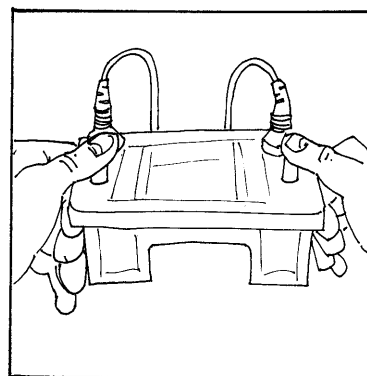- Gently lift the gel tray from the unit. **Always wear gloves, eye protection and protective clothing** if buffer and/or gel contains Ethidium Bromide. Ethidium Bromide is a powerful mutagen, gloves, eye protection and protective clothing should always be worn when handling the gel or buffer solutions. See Material Data Safety Sheets.

- View separated fragments under UV light, using proper protection for eyes and skin (see manufacturer's instructions).

3.4 Maximum Well/Comb Volumes

NOTE: To calculate sample well volume expressed in millimeters (mm) of height, divide maximum volume by tooth depth.

| MBU Comb for MGU-402T # of wells | Tooth Depth mm | Overall length of comb mm | Spacing between teeth mm | Tooth width/mm | 1.0mm thickness volume per tooth microliters (ul) | 1.5mm thickness volume per tooth microliters (ul) | 2.0mm thickness volume per tooth microliters (ul) | 3.0mm thickness volume per tooth microliters (ul) |
|---|----------------|---------------------------|--------------------------|----------------|---|---|---|---|
| 5 | 12.7 | 87.28 | 1.58 | 16.2 | 205.74 | 308.61 | 411.48 | 617.22 |
| 12 | 12.7 | 87.28 | 1.58 | 5.83 | 74.04 | 111.06 | 148.08 | 222.12 |
| 18 | 12.7 | 87.28 | 1.58 | 3.37 | 42.79 | 64.19 | 85.58 | 128.37 |
| SGC14 Comb for MGU-502T & MGU-602T # of wells | Tooth Depth mm | Overall length of comb mm | Spacing between teeth mm | Tooth width/mm | 1.0mm thickness volume per tooth microliters (ul) | 1.5mm thickness volume per tooth microliters (ul) | 2.0mm thickness volume per tooth microliters (ul) | 3.0mm thickness volume per tooth microliters (ul) |
| 1 | 12.7 | 120.64 | 3 | 120 | 1524 | 2286 | 3048 | 4672 |
| 3E | 12.7 | 120.64 | 3 | 38.1 | 483.87 | 725.81 | 967.74 | 1451.61 |
| 5 | 12.7 | 120.64 | 3 | 21.6 | 274 | 411.48 | 548 | 822 |
| 10 | 12.7 | 120.64 | 3 | 9.2 | 117 | 175.26 | 232 | 350 |
| 12 | 12.7 | 120.64 | 2.4 | 7.8 | 99 | 148.59 | 198 | 297 |
| 13 | 12.7 | 120.64 | 2.4 | 6.6 | 83.82 | 125.73 | 167.64 | 251.46 |
| 16 | 12.7 | 120.64 | 2.4 | 5.3 | 67.3 | 100.9 | 134.6 | 201 |
| 20 | 12.7 | 120.64 | 1.6 | 4.5 | 57 | 85.7 | 114 | 171 |
| 26 | 12.7 | 120.64 | 1.6 | 3.1 | 39.3 | 59 | 78 | 118 |
| 27 | 12.7 | 120.64 | 1.6 | 2.93 | 37.21 | 55.82 | 74.42 | 111.63 |

SECTION 4

Running Conditions

4.1 Recommended Power

As a rule, optimal resolution of larger molecules is achieved during longer runs at lower voltages, whereas smaller molecules require shorter runs at higher voltages. Applied voltage gradients can therefore be anywhere in the range of 1-10 VDC/cm of gel. Using the standard buffer systems listed below, most runs will use 5 VDC/cm as a general rule.

The usual run time will vary for the voltage chosen but should range from 15 to 60 minutes. Nucleic acid migration is monitored by the progress of marker dyes. Constant power is not a necessity, but it produces uniform heat throughout the run, therefore minimizing band diffusion. Be sure the polarity is correct i.e. that the DNA is loaded near the cathode (black electrode) to run toward the anode (red terminal).

Agarose gels may be stored for several days at 4°C wrapped in plastic wrap. Seakem Agarose (FMC) is used (normally) for preparative and analytical gels. Other types of agarose can be used for special purposes.

4.2 Recommended Buffers*

| Type* | Concentrated Stock/liter | Final Concentration |
|------------------------------------|---|--|
| TAE (Tris-acetate) | 50X - 242 gm Tris base 57.1ml glacial acetic acid 100 ml 0.5M EDTA (pH 8.0) | 1X - 0.04M Tris-acetate 0.001M EDTA |
| TBE (Tris-borate) | 5X - 54 gm Tris base 27.5 gm boric acid 20 ml 0.5M EDTA (pH 8.0) | 0.5X - 0.045M Tris-borate 0.001M EDTA |
| <u>10X - Loading Buffer (DNA)*</u> | | |
| 0.25% Bromophenol blue | | |
| 0.25% Xylene cyanol | | |
| 20% Ficoll Type 400 | | |
| 0.1M EDTA, pH 8.0 | | |

Ethidium Bromide Staining

EtBr can be premixed with buffers and agarose for use during electrophoresis. Add to agarose only after temperature has fallen below 55°C. Gels can also be stained after electrophoresis in a soaking tray. Use EtBr at a final concentration of 0.1 µg/ml from a 1mg/ml stock solution. Ethidium Bromide is a powerful mutagen. ALWAYS wear gloves, eye protection and protective clothing. Dispose of solutions in accordance with the safety regulations of your institution.

4.3 References

1. Hames, B.D., Rickwood, D. (ed.) (1990). *Gel Electrophoresis of Nucleic Acids. A Practical Approach*. 2nd edn. IRL Press, Oxford. Ch. 2.
2. Sambrook, J., Fritsch, E.F., Maniatis, T. (1989). *Molecular Cloning. A Laboratory Manual*. 2nd edn. Cold Spring Harbor Laboratory Press, Cold Spring Harbor, New York. Ch 6.
3. Ausubel, F.M., Brent, R., Kingston, R.E., Moore, D.D., Seidman, J.G., Smith, J.A., Struhl, K. (ed) (1993). *Current Protocols in Molecular Biology*. Vol. 1, Greene Publishing Associates, Inc. and John Wiley & Sons, Inc., Ch. 2.

SECTION 5 Maintenance of Equipment

5.1 Care and Handling

The plastic components of the Midi-Horizontal units are fabricated from acrylic and polycarbonate. Electrodes and connectors are made from pure platinum, stainless steel, and chrome plated brass. As with any laboratory instrument, adequate care ensures consistent and reliable performance.

After each use, rinse buffer chamber, gel tray and combs with de-ionized water. Wipe dry with a soft cloth or paper towel, or allow to air dry. Whenever necessary, all components may be washed gently with water and a non-abrasive detergent, and rinsed and dried as above. *Never* use abrasive cleaners, glass cleaning sprays or scouring pads to clean the components, as these will damage the unit and components.

Additional precautions:

- Do not autoclave or dry-heat sterilize the apparatus or components.
- Do not expose the apparatus or components to phenol, acetone, benzene, halogenated hydrocarbon solvents or alcohols.
- Avoid prolonged exposure of the apparatus or components to UV light.
- Do NOT treat with diethylpyrocarbonate (DEPC)-treated water for extended periods at 37°C. A brief rinse with DEPC-water is sufficient after a thorough wash.

5.2 Maintenance

The following inspection and maintenance procedures will help maintain the safety and reliable performance of the Midi-Horizontal systems. Replacement parts can be ordered by calling

1-858-755-4959 or by contacting your local distributor.

- Banana plugs and power cords should be inspected regularly. If the banana plugs become loose or do not feel friction tight replace the plugs or power cords.
- Should power cord assemblies (connectors, wire or shrouds) show any signs of wear or damage (e.g. cracks, nicks, abrasions, or melted insulation), replace them immediately.
- The platinum wire is secured to the banana jack by compression between a stainless washer and the jack nut. The nut/washer interface should be tight and free of corrosion.

SECTION 6

Horizontal Midi-Gel Kits and Accessories

| Cat.# | Item |
|-------------------|---|
| GCMGU-402T | Horizontal Midi-Gel Productivity Package CE , with tapered baffles. Gel bed dimensions are 11cm x 11cm. Includes two combs, two UVT gel tray, external gel caster and safety cover. Please specify comb cat. # from list below. |
| GCMGU-502T | Double-Wide Midi-Gel Productivity Package CE , with tapered baffles. Gel bed dimensions are 14cm x 10cm. Includes two combs, two UVT gel tray, external gel caster and safety cover. Please specify comb cat. # from list below. |
| GCMGU-602T | Horizontal Midi-Gel Productivity Package CE , with tapered baffles. Gel bed dimensions are 14cm x 14cm. Includes two combs, two UVT gel tray, external gel caster and safety cover. Please specify comb cat. # from list below. |

Color Coded Midi-Gel Option:

Available for the MGU-402T, MGU-502T and MGU-602T models. To specify color desired, please add the following color code to the end of the catalog number: -Y (Yellow), -P (Purple), -B (Blue), -PK (Pink), or -G (Green). For example, to specify a green MGU-402T place order for catalog # MGU-402T-G.

MIDI-GEL ACCESSORIES

MGU-402T combs (Available in Polycarbonate, or Teflon-Coated Aluminum)**

| Cat.# | Comb Dimensions |
|-------|-----------------|
|-------|-----------------|

| | |
|----------|-------------------|
| MBU-1005 | 1mm x 5 well |
| MBU-1010 | 1mm x 10 well** |
| MBU-1012 | 1mm x 12 well |
| MBU-1018 | 1mm x 18 well |
| MBU-1505 | 1.5mm x 5 well |
| MBU-1510 | 1.5mm x 10 well** |
| MBU-1512 | 1.5mm x 12 well |
| MBU-1518 | 1.5mm x 18 well |

** To specify Teflon combs add "-TF" to end of cat. #.

MGU-502T and MGU-602T combs (Available in Polycarbonate, or Teflon-Coated Aluminum)**

| Cat.# | Comb Dimensions |
|-------|-----------------|
|-------|-----------------|

| | |
|-------------|--------------------------------|
| SGC14-1001 | 1mm x 1 well |
| SGC14-1003E | 1mm x 3 equal wells |
| SGC14-1003M | 1mm x 2 well + middle marker |
| SGC14-1005 | 1mm x 5 well |
| SGC14-1008 | 1mm x 8 well |
| SGC14-1010 | 1mm x 10 well |
| SGC14-1012 | 1mm x 12 well |
| SGC14-1013 | 1mm x 13 well* |
| SGC14-1016 | 1mm x 16 well |
| SGC14-1020 | 1mm x 20 well |
| SGC14-1026 | 1mm x 26 well |
| SGC14-1027 | 1mm x 27 well* |
| SGC14-1501 | 1.5mm x 1 well |
| SGC14-1503E | 1.5mm x 3 equal wells |
| SGC14-1503M | 1.5mm x 2 well + middle marker |
| SGC14-1505 | 1.5mm x 5 well |
| SGC14-1508 | 1.5mm x 8 well |
| SGC14-1510 | 1.5mm x 10 well |
| SGC14-1512 | 1.5mm x 12 well |
| SGC14-1513 | 1.5mm x 13 well* |
| SGC14-1516 | 1.5mm x 16 well |
| SGC14-1520 | 1.5mm x 20 well |
| SGC14-1526 | 1.5mm x 26 well |
| SGC14-1527 | 1.5mm x 27 well |

| Cat.# | Comb Dimensions |
|-------|-----------------|
|-------|-----------------|

| | |
|-------------|------------------------------|
| SGC14-2001 | 2mm x 1 well |
| SGC14-2003E | 2mm x 3 equal wells |
| SGC14-2003M | 2mm x 2 well + middle marker |
| SGC14-2005 | 2mm x 5 well |
| SGC14-2008 | 2mm x 8 well |
| SGC14-2010 | 2mm x 10 well |
| SGC14-2012 | 2mm x 12 well |
| SGC14-2013 | 2mm x 13 well* |
| SGC14-2016 | 2mm x 16 well |
| SGC14-2020 | 2mm x 20 well |
| SGC14-2026 | 2mm x 26 well |
| SGC14-2027 | 2mm x 27 well* |
| SGC14-3001 | 3mm x 1 well |
| SGC14-3003E | 3mm x 3 equal wells |
| SGC14-3003M | 3mm x 2 well + middle marker |
| SGC14-3005 | 3mm x 5 well |
| SGC14-3008 | 3mm x 8 well |
| SGC14-3010 | 3mm x 10 well |
| SGC14-3012 | 3mm x 12 well |
| SGC14-3013 | 3mm x 13 well* |
| SGC14-3016 | 3mm x 16 well |
| SGC14-3020 | 3mm x 20 well |
| SGC14-3026 | 3mm x 26 well |
| SGC14-3027 | 3mm x 27 well |

*Designed for multi-channel pipettor loading.

** To specify Teflon combs add "-TF" to end of cat. #.

MIDI-GEL ACCESSORIES-CONTINUED

Midi Gel Trays (UV Transparent)

| Cat. # | Item |
|--------|------|
|--------|------|

| | |
|---------|--|
| MGT-400 | Gel Tray for the MGU-402T, 2 slots. Pathlength of 4cm/slot |
| MGT-500 | Gel Tray for the MGU-502T, 2 slots. Pathlength of 3.6cm/slot |
| MGT-600 | Gel Tray for the MGU-602T, 4 slots. Pathlength of 3.0cm/slot |

Replacement Baffles

| Cat. # | Item |
|--------------|------------------------------|
| BAFFLES-402T | Tapered Baffles for MGU-402T |
| BAFFLES-502T | Tapered Baffles for MGU-502T |
| BAFFLES-602T | Tapered Baffles for MGU-602T |

Horizontal Gel Casters

| Cat. # | Item |
|-----------|--|
| GCT10-111 | Gel Casting Tray for 10.5cm (w) x 11cm (l), single place |
| GCT10-112 | Gel Casting Tray for 10.5cm (w) x 11cm (l), two place |
| GCT14-101 | Gel Casting Tray for 14cm (w) x 10cm (l), single place |
| GCT14-102 | Gel Casting Tray for 14cm (w) x 10cm (l), two place |
| GCT14-161 | Gel Casting Tray for 14cm (w) x 16cm (l), single place |
| GCT14-162 | Gel Casting Tray for 14cm (w) x 16cm (l), two place |

Gel Tape

| Cat. # | Item |
|----------|--------------------------|
| GT-72-10 | Gel tape, 72 yard x 1.0" |
| GT-72-15 | Gel tape, 72 yard x 1.5" |

Power Supply

| Cat. # | Item |
|--------------|---|
| EPS-200-II | Mini Power Supply, 24 hour timer, CV or CC, 10-200V, 110V/60Hz, current range:4-2000mA, 200 watts |
| EPS-200-IIV | Mini Power Supply, 24 hour timer, CV or CC, 10-200V, 220V/50Hz, current range:4-2000mA, 200 watts |
| EPS-300-II | Mini Power Supply, 24 hour timer, CV or CC, 10-300V, 110V/60Hz, current range:4-500mA, 90 watts |
| EPS-300-IIV | Mini Power Supply, 24 hour timer, CV or CC, 10-300V, 220V/50Hz, current range:4-500mA, 90 watts |
| EPS-3000-III | Constant Power Supply, 0-3000 Volt, 110V/60Hz, Series II |
| EPS-3000-IIV | Constant Power Supply, 0-3000 Volt, 220V/50Hz, Series II |

Replacement Safety Covers with attached power leads

| Cat. # | Item |
|-------------|--|
| MGU-402T-SC | Replacement safety cover with power leads for MGU-402T |
| MGU-502T-SC | Replacement safety cover with power leads for MGU-502T |
| MGU-602T-SC | Replacement safety cover with power leads for MGU-602T |

NOTES

NOTES

CONTACT US

Telephone:
Local or International
858-755-4959
Toll Free: 858-755-0733

Fax: 858-755-0733

Check out our new and improved online ordering!

www.cbsscientific.com

E-mail address:
sales@cbssci.com

Mailing address:
C.B.S. Scientific Company
P.O. Box 856
Del Mar, CA 92014

Shipping address:
C.B.S. Scientific Company
10805 Vista Sorrento Pkwy
Suite 100
San Diego, CA 92121

Credit Card Options:
Visa/Mastercard
Discover/American Express