
# INSTRUCTION MANUAL

## **Pore Gradient Lipoprotein Electrophoresis System (LPE) LPE-4003**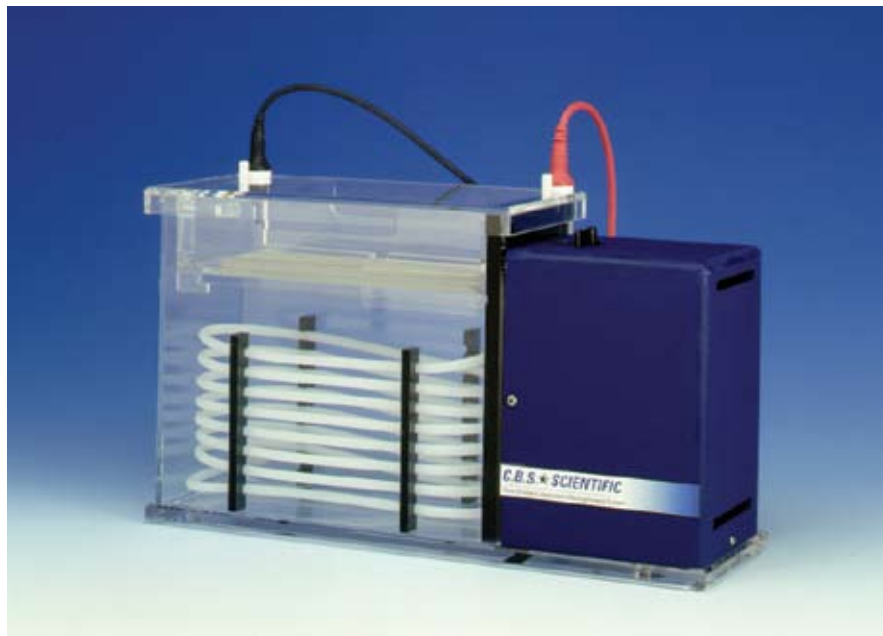

# TABLE OF CONTENTS

	Page
<b>Important User Information</b>	3-4
<b>Section 1    General Information</b>	
1.1    Introduction and Features	5
1.2    Specifications	5
1.3    Safety	5
<b>Section 2    Description of parts</b>	
2.1    Unpacking	6
2.2    Components/Assembly	6
<b>Section 3    Instructions for Use</b>	
3.1    PGGE Unit Preparation	7
3.2    Buffer Preparation	7
3.3    Gel Preparation when using pre-cast gels	7
3.4    Running Standard PAGE gels	7
3.5    Sample Loading and Electrophoresis	7
3.6    Gel Removal	7
<b>Section 4    Running Conditions</b>	
4.1    Recommended Power	7
4.2    References	7
<b>Section 5    Maintenance of Equipment</b>	
5.1    Care and Handling	8
5.2    Maintenance	8
<b>Section 6    LPE Replacement parts list</b>	9

## IMPORTANT USER INFORMATION

This Instruction Manual will explain how to use this product safely and effectively. Please read and carefully follow the instruction manual in its entirety.

The triangle/exclamation mark symbol alerts the user of the product to important operational, maintenance, and/or warranty requirements.

The triangle/lighting bolt symbol alerts the user of the product to potentially hazardous electrical exposure.

Failure to adhere to the instructions could result in personal and/or laboratory hazards, as well as invalidate any warranty. Always turn off the DC power source prior to disconnecting power cords from the product. Disconnect power cords from the power source first and then from the product. For maximum safety, always operate this system in an isolated, low traffic area, not accessible to unauthorized personnel. Never operate damaged or leaking equipment.

## WARRANTY AND LIABILITY

This product was produced utilizing the highest practical standards of materials, workmanship, and design. C.B.S. SCIENTIFIC warrants that the product has been tested and will meet or exceed published specifications. This warranty is valid only if the product has been operated and maintained according to the instructions provided.

C.B.S. SCIENTIFIC warrants this product to be free from defects in materials and workmanship under normal service for one year from date of shipment. If the product proves defective during this period, SCIENTIFIC will repair or replace it at our option, free of charge, if returned to us postage prepaid. This warranty does not cover: damage in transit, damage caused by carelessness, misuse or neglect, normal wear through frequent use, damage caused by solvent corrosion, damage caused by improper handling or user alteration, nor unsatisfactory performance as a result of conditions beyond our control. C.B.S. SCIENTIFIC shall in no event be liable for incidental nor consequential damages, including without limitation, lost profits, loss of income, loss of business opportunities, loss of use and other related damages, however caused, nor any damage arising from the incorrect use of the product.

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## SECTION 1 General Information

### 1.1 Introduction

The PGGE, Pore Gradient Lipoprotein Electrophoresis System, Cat. # LPE-4003, is designed to run 1-4 nondenaturing gradient gels with dimensions of 82 x 82 x 5mm (Alamo Gels). The primary application is for the resolution of Lipoprotein sub-classes (LDL, HDL) based on particle size using non-denaturing pore gradient gel electrophoresis (GGE). Other applications include Standard PAGE and Short DNA Sequencing runs. The PGGE system features; 4-gel capacity, temperature control, buffer recycling and inter-locking safety cover.

**Table 1  
Features of PGGE System**

Model #		Glass Plate Dimensions (mm) (w x h)	Precast Gels available	Internal cooling chamber	Manual Gel Casting	Buffer Cycling
LPE-4003	4 gel capacity	82 x 82 x 5	Yes	Yes	Yes	Yes

### 1.2 Specifications

**Constructions:**

**Buffer chamber, safety cover**

**Electrodes**

**Power cords**

**Safety Certification**

Acrylic, Kydex

Platinum wire .012" diameter

FR Urethane rated 7500VDC, 200mA, 65°C

EN61010-1-1993 (IEC1010-1)

**Table 2  
Specifications**

	LPE-4003
Shipping weight	12 lbs.
Overall Size (l)x(w)x(h) cm	39 x 12 x 28
Recommended buffer volume	3.5 L
Distance between electrodes -cm	16
Voltage limit	250VDC
Mains	115V/60Hz

### 1.3 Safety

Power to the PGGE unit is to be supplied by an external DC voltage power supply that must be ground isolated so that the DC voltage output floats with respect to ground. For any power supply used, the maximum specified operating parameters for the units are:

**Maximum Limits**

250VDC

30watts power

125mA current

55°C ambient temperature

Current to the unit, provided from the external power supply, must enter the unit through the lid assembly, providing a safety interlock to the user. Current to the unit is broken when the lid is removed. **Do not attempt to use the unit without the safety lid, and always turn the power supply off before removing the lid, or when working with the unit in any way. Follow safety precautions specified by the power supply manufacturer.**

## SECTION 2


### Description of Parts

#### 2.1 Unpacking

Please verify that your unit comes complete with the following components:

- Base unit with lower buffer chamber, cooling coil with external connections and recycling pump
- Upper reservoir with 2 grey silicone gaskets
- Safety Cover & power leads
- 3 gel slot plugs
- Mains-power cord (not shown)

#### 2.2 Components/Assembly

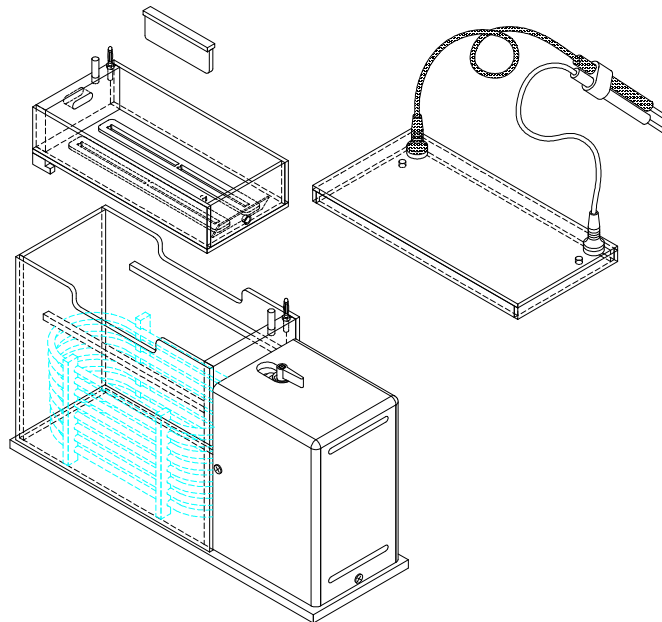## SECTION 3

### Instructions for Use

#### 3.1 PGGE Unit Preparation

Place the unit on a level work surface in an authorized work area. Attach the black (Mains) power cord to the unit and a mains outlet. Check to be sure the toggle switch is in the "0" position.

#### 3.2 Buffer Preparation

Prepare the PGGE buffer as in Ref. 1 (90mM Tris, 9mM borate, 2.5mM EDTA, pH 8.3). The buffer should be pre-chilled to 5°C (cold room). Add 3.5L of chilled buffer to the lower reservoir. Connect the inlet and outlet tubes to the refrigerated water bath (chiller). For optimal running conditions, set the waterbath temperature control to 7°C and begin circulation.

#### 3.3 Gel Preparation when using pre-cast gels

Remove the pre-cast gel cassettes from protective wrapping (always wear gloves!). Rinse in d.d.H<sub>2</sub>O. Insert gel cassettes into upper reservoir such that 1cm remains above the grey gasket. Turn toggle switch to "I" position to begin buffer cycling and filling upper reservoir to correct level.

#### 3.4 Running Standard PAGE gels.

The LPE-4003 can be used for other types of gels other than pre-cast non-denaturing acrylamide gradients. Accessories for manually casting PAGE gels are available for the LPE-4003. Please contact our customer service department for ordering information.

#### 3.5 Sample Loading and Electrophoresis

Prepare samples as per ref. 1,2,3. Turn buffer cycling off, insert well former into gel cassettes and load samples with controls. Resume buffer cycling, attach safety cover, connect power cords to DC power supply and begin electrophoresis.

#### 3.6 Gel Removal

After completion of electrophoresis, turn off power supply and disconnect power cords. Remove the upper reservoir. Remove gel cassettes from grey gasket by pushing from top until cassette ejects from gasket.

## SECTION 4

### Running Conditions

#### 4.1 Recommended Power

Precise electrophoresis conditions will vary according to the number and type of gels used, buffer conditions employed, power input, and the general goal of the experiment. Refer to the reference section for in depth discussions on practical; and theoretical approaches to Lipoprotein electrophoresis.

#### 4.2 References

Rainwater, David L., et.al. Production of polyacrylamide gradient gels for the electrophoretic resolution of lipoproteins. *Journal of Lipid Research*, **Volume 33**, 1992.

Rainwater, David L., Electrophoretic Separation of LDL and HDL Subclasses, *Methods in Molecular Biology*, **Volume 110**: Lipoprotein Protocols.

Rainwater, David L., et.al. Characterization of a composite gradient gel for the electrophoretic separation of lipoproteins, *Journal of Lipid Research*, **Volume 38**, 199.

## SECTION 5

### Maintenance of Equipment

#### 5.1 Care and Handling

The components of the PGGE units are fabricated from acrylic, Kydex, silicone rubber and polyethylene. Electrodes and connectors are made from pure platinum, stainless steel, and chrome plated brass. As with any laboratory instrument, adequate care ensures consistent and reliable performance.

After each use, rinse buffer chamber and accessories with de-ionized water. Wipe dry with a soft cloth or paper towel, or allow to air dry. Whenever necessary, all components may be washed gently with water and a non-abrasive detergent, and rinsed and dried as above. *Never* use abrasive cleaners, glass cleaning sprays or scouring pads to clean the components, as these will damage the unit and components.

Additional precautions:

- Do not autoclave or dry-heat sterilize the apparatus or components.
- Do not expose the apparatus or components to phenol, acetone, benzene, halogenated hydrocarbon solvents or alcohols.
- Avoid prolonged exposure of the apparatus or components to UV light.
- Do NOT treat with diethylpyrocarbonate (DEPC)-treated water for extended periods at 37°C. A brief rinse with DEPC-water is sufficient after a thorough wash.

#### 5.2 Maintenance


The following inspection and maintenance procedures will help maintain the safety and reliable performance of the PGGE systems. Replacement parts can be ordered by calling 1-800-243-4959 or by contacting your local distributor.

- Banana plugs and power cords should be inspected regularly. If the banana plugs become loose or do not feel friction tight replace the plugs or power cords.
- Should power cord assemblies (connectors, wire or shrouds) show any signs of wear or damage (e.g. cracks, nicks, abrasions, or melted insulation), replace them immediately.
- The platinum wire is secured to the banana jack by compression between a stainless washer and the jack nut. The nut/washer interface should be tight and free of corrosion.

## SECTION 6

### LPE-4002 REPLACEMENT PARTS LIST

Part Number	Qty/unit	Description
LPE-4002-001	2	Grommet 2x80
LPE-4002-002	3	Plug for 2X80 grommet
LPE-4002-003	1	Pump
LPE-4002-004	1	Safety Cover W/Power Leads
LPE-4002-005	1	Tank, upper
LPE-4002-006	1	Tank, lower
LPE-4002-007	1	Pump cover w/AC switch/outlet
LPE-4002-008	1	1/8" NPT Nipple
LPE-4002-009	1	Hose, 5"
LPE-4002-010	1	Plastic ball valve
LPE-4002-011	1	Grommet, pump intake
LPE-4002-012	1	polytube
LPE-4002-013	1	RA fitting
LPE-4002-014	1	O-Ring

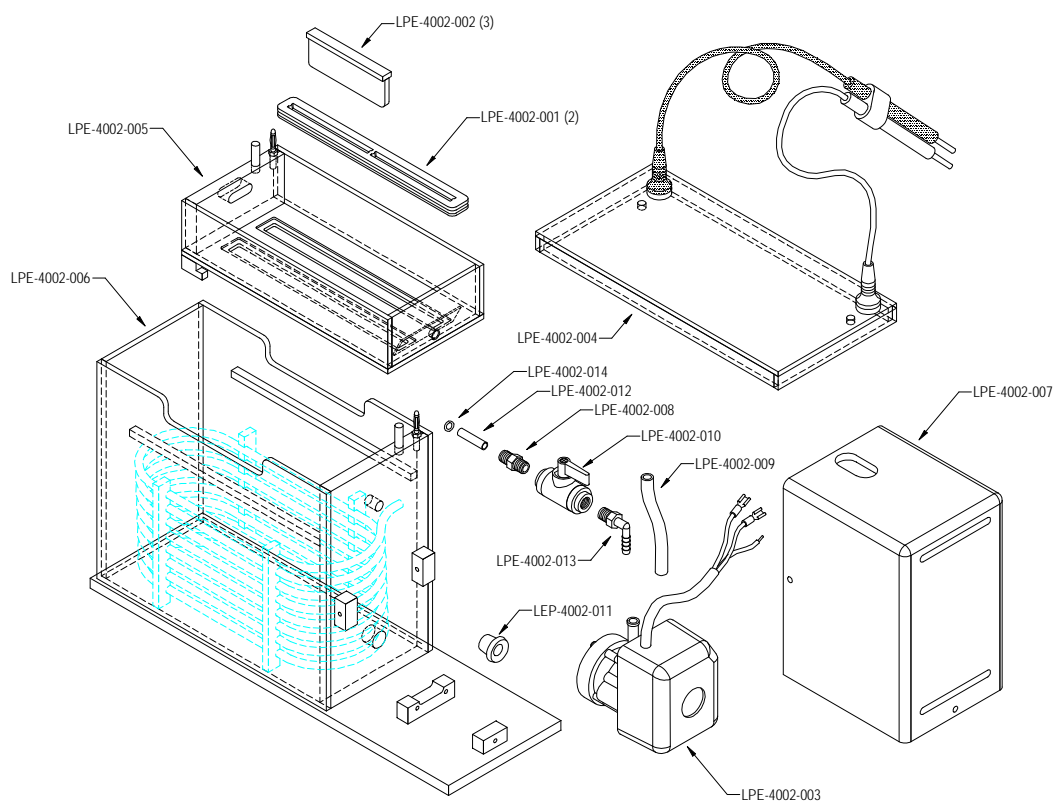#### POWER SUPPLY

Cat. #	Item
EPS-250	Mini Power Supply, 20-250 Volts, 300mA @ 150-250V or 500mA below 150V
EPS-250V	Mini Power Supply, 20-250 Volts, 300mA @ 150-250V or 500mA below 150V, 220V
EPS-250-II	Mini Power Supply, 10-250 Volts, 10-500mA with timer
EPS-250-IIV	Mini Power Supply, 10-250 Volts, 10-500mA with timer, 220V

---

---

## NOTES

---

---

---

---

## NOTES

---

---

# CONTACT INFORMATION

**Telephone:**  
Local or International  
858-755-4959  
Toll Free: 800-243-4959

**Fax:** 858-755-0733

**Online ordering:**  
[www.cbsscientific.com](http://www.cbsscientific.com)

**E-mail address:**  
[sales@cbssci.com](mailto:sales@cbssci.com)

**Mailing address:**  
C.B.S. Scientific Company  
P.O. Box 856  
Del Mar, CA 92014

**Shipping address:**  
10805 Vista Sorrento Parkway  
Ste 100  
San Diego, CA 92121

**Credit Card Options:**  
Visa/Mastercard  
Discover/American Express