
# INSTRUCTION MANUAL

## Horizontal Mini-Gel Systems

**MGU-102T**

**MGU-202T**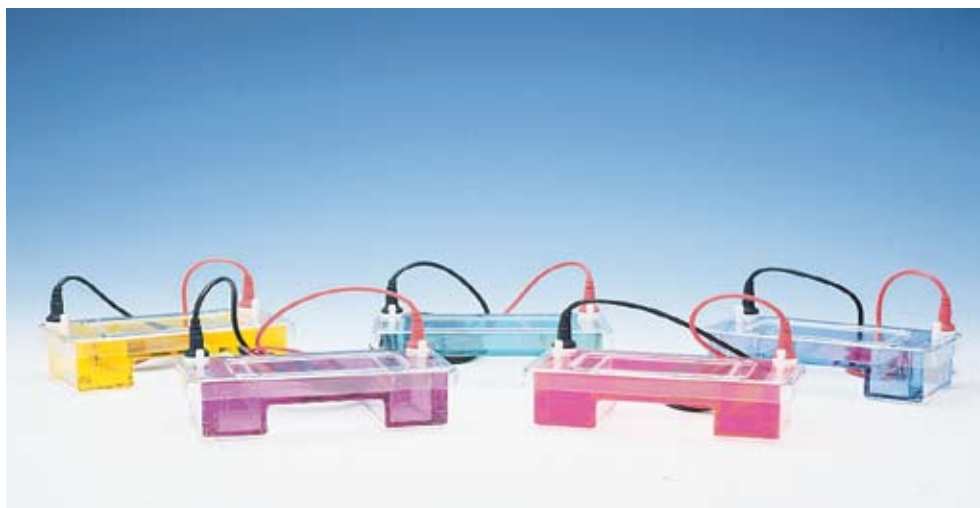

# TABLE OF CONTENTS

	Page
<b>Important User Information</b>	<b>3-4</b>
<b>Section 1    General Information</b>	
1.1    Introduction	5
1.2    Specifications	5
1.3    Safety	5
<b>Section 2    Description of parts</b>	
2.1    Unpacking	6
2.2    Components/Assembly	6
<b>Section 3    Instructions for Use</b>	
3.1    Mini-Gel Unit Preparation and Gel Casting	7-8
3.2    Running the Gel	9
3.3    Removing the Gel	9
3.4    Maximum well/Comb volumes	10
<b>Section 4    Running Conditions</b>	
4.1    Recommended Power	10
4.2    Recommended Buffers	10-11
<b>Section 5    Maintenance of Equipment</b>	
5.1    Care and Handling	11
5.2    Maintenance	11
<b>Section 6    Horizontal Mini-Gel Kits and Accessories</b>	<b>12</b>

## IMPORTANT USER INFORMATION

This Instruction Manual will explain how to use this product safely and effectively. Please read and carefully follow the instruction manual in its entirety.

The triangle/exclamation mark symbol alerts the user of the product to important operational, maintenance, and/or warranty requirements.

The triangle/lightning bolt symbol alerts the user of the product to potentially hazardous electrical exposure.

Failure to adhere to the instructions could result in personal and/or laboratory hazards, as well as invalidate any warranty. Always turn off the DC power source prior to disconnecting power cords from the product. Disconnect power cords from the power source first and then from the product. For maximum safety, always operate this system in an isolated, low traffic area, not accessible to unauthorized personnel. Never operate damaged or leaking equipment.

## WARRANTY AND LIABILITY

This product was produced utilizing the highest practical standards of materials, workmanship, and design. C.B.S. Scientific warrants that the product has been tested and will meet or exceed published specifications. This warranty is valid only if the product has been operated and maintained according to the instructions provided.

C.B.S. Scientific warrants this product to be free from defects in materials and workmanship under normal service for one year from date of shipment. If the product proves defective during this period, C.B.S. Scientific will repair or replace it at our option, free of charge, if returned to us postage prepaid. This warranty does not cover: damage in transit, damage caused by carelessness, misuse or neglect, normal wear through frequent use, damage caused by solvent corrosion, damage caused by improper handling or user alteration, nor unsatisfactory performance as a result of conditions beyond our control. C.B.S. Scientific shall in no event be liable for incidental nor consequential damages, including without limitation, lost profits, loss of income, loss of business opportunities, loss of use and other related damages, however caused, nor any damage arising from the incorrect use of the product.

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### General Information

#### 1.1 Introduction

The Horizontal Mini-Gel systems, MGU-102T and MGU-202T, provide a compact format to rapidly separate and identify DNA and RNA fragments, PCR products and synthetic oligonucleotides, while using minimal quantities of expensive reagents. The tapered baffle system allows gel casting directly onto the unit or removable trays without having to seal with agarose plugs or tape. The gel bed and gel tray are UV transparent. Combs are offered in several thicknesses and formats to accommodate a wide variety of experimental applications.

#### 1.2 Specifications

##### Constructions

Buffer chamber, lid	Acrylic
Electrodes	Platinum wire .012" diameter
Power cords	FR Urethane, rated 7500V, 200mA, 65° C
Tapered baffles	Acrylic
Combs	Polycarbonate, Anodized Aluminum or Teflon
Adjustable comb backing	Acrylic
UVT Gel Tray (optional)	U-V transparent acrylic
Safety Certification	EN61010-1-1993 (IEC1010-1)

	MGU-102T	MGU-202T
Shipping weight	3 lb.	3 lb.
Overall Size (l)x(w)x(h)	18.2cm x 8cm x 9cm	20cm x 9.7cm x 10cm
Gel Size	5.5cm x 8.5cm	7.5cm x 10cm
Recommended buffer volume:		
with gel tray	150 ml	250 ml
without gel tray	140 ml	235 ml
Distance between electrodes	14.5 cm	21 cm
Voltage limit	250VDC	250VDC

#### 1.3 Safety

Power to the MGU-102T or MGU-202T units is to be supplied by an external DC voltage power supply that must be ground isolated so that the DC voltage output floats with respect to ground. For any power supply used, the maximum specified operating parameters for the units are:

##### Maximum Limits

250VDC voltage  
15 watts power  
60 mA current  
50°C ambient temperature

Current to the unit, provided from the external power supply, must enter the unit through the lid assembly, providing a safety interlock to the user. Current to the unit is broken when the lid is removed. **Do not attempt to use the unit without the safety lid. Always turn the power supply off before removing the lid, or when working with the unit in any way. Follow safety precautions specified by the power supply manufacturer.**

## SECTION 2


## Description of Parts

### 2.1 Unpacking

Please verify that your MGU-102T or MGU-202T unit comes complete with the following components:

- Mini-Horizontal electrophoresis unit
- Tapered baffles (2)
- Safety cover with attached DC power cords
- Adjustable height comb (1 each of your choice)

### 2.2 Components/Assembly

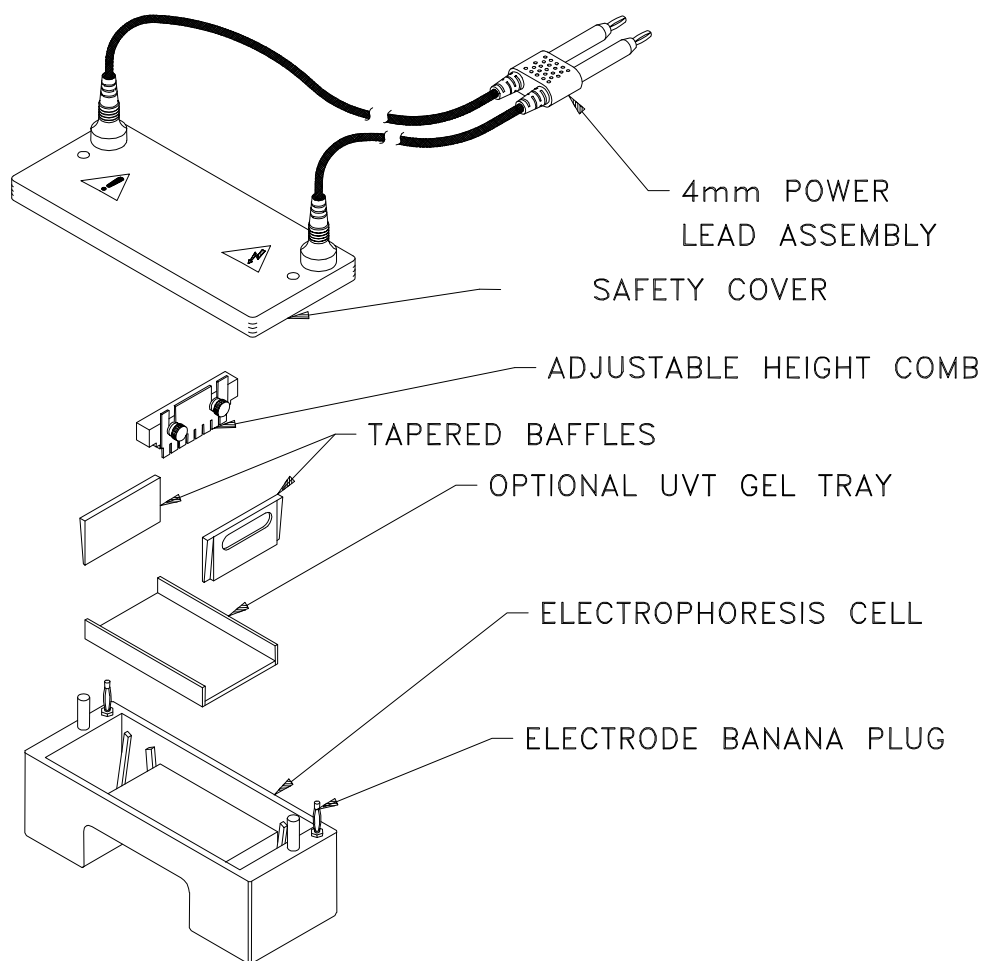## SECTION 3

### Instructions for Use


#### 3.1 Mini-Gel Unit Preparation and Gel Casting

Place Mini-Gel unit on a level work surface in an authorized work area. Place the adjustable height comb assembly over the gel bed or gel tray. Loosen the white screws holding the comb to the backing and adjust comb depth (generally 2mm). Gently tighten the adjusting screws and set the comb aside. **See Table 1 for comb volumes.**


The Mini-Gel unit provides the user with three options for gel casting: (a) **gel casting directly into unit**, (b) **gel casting directly into unit using gel tray**, or (c) **gel casting in gel tray outside of unit**.

##### A. Gel Casting Directly into Unit.

1. Slide the tapered baffles snugly into the grooves next to the bed, handles to the outside. The tapered baffles will rest friction tight against the gel bed. Do NOT press the baffles too tightly into the stops.

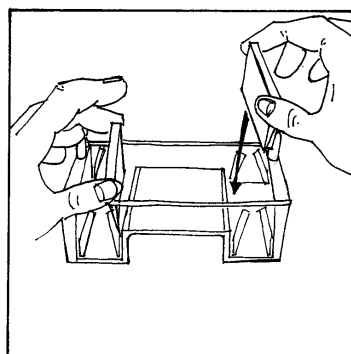
2. Position the gel comb, remembering, current travels from negative (black cathode) towards positive (red anode). Pour the agarose solution onto the gel bed. Allow the agarose slab to cool uncovered until it turns opaque.
3. Slide the baffles out of their slots and store in the slots provided at either end of the unit near both terminals.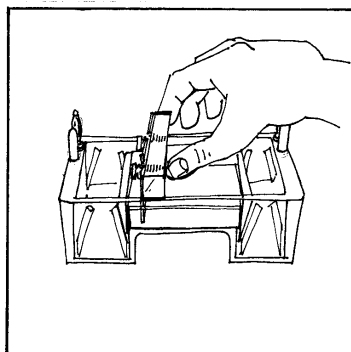


##### B. Casting Directly into Unit Using Gel Tray.

1. Insert gel tray in unit. Simultaneously slide the tapered baffles snugly into the grooves next to the bed, handles to the outside at the same time. The tapered baffles will rest friction tight against the gel tray. Do NOT press the baffles too tightly into the stops.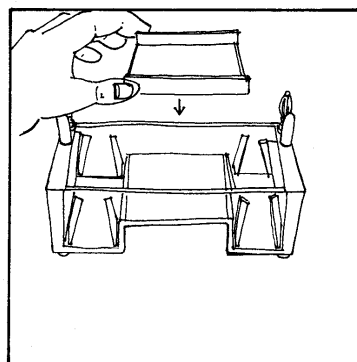



2. Position the gel comb, remembering, current travels from negative (black cathode) towards positive (red anode). Pour the agarose solution onto the gel bed. Allow the agarose slab to cool uncovered until it turns opaque.
3. Slide the baffles out of their slots and store in the slots provided at either end of the unit neat both terminals.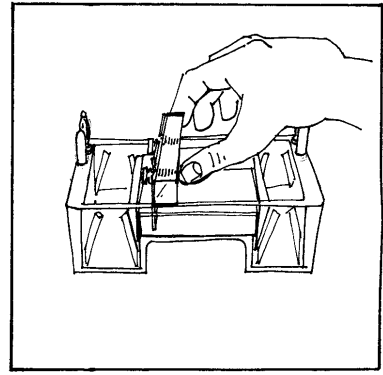


**C. Gel Casting into Gel Tray Outside of Unit.**

1. Tape both ends of gel tray with tape.
2. Position the gel comb in the gel tray at one end. Pour the agarose solution onto the gel bed. Allow the agarose slab to cool uncovered until it turns opaque.  
NOTE: (Optional storage) Pre-cast gels can be stored for up to one month if wrapped in plastic wrap and stored at 4°C. Remove combs prior to storing.

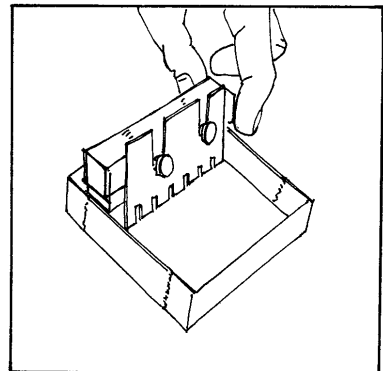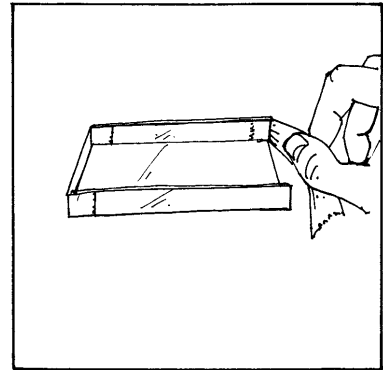
3. Remove comb and gel tape at each end. Position the tray in the unit, remembering, current travels from negative (black cathode) towards positive (red anode).
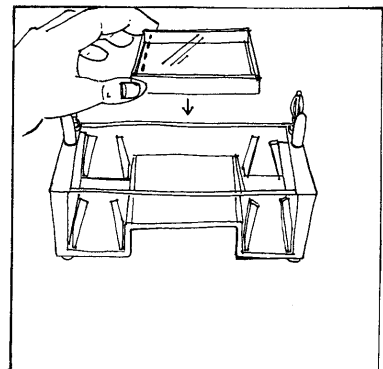


### 3.2 Running the gel

1. Add enough buffer to fill both reservoirs and overflow the surface of the gel to a depth of 2-3mm. Gently remove the comb. Flush out any air bubbles in the wells. Load the samples into sample wells. **NOTE: DO NOT FORGET TO LOAD DNA SIZE STANDARD.** Place the safety cover on the unit.

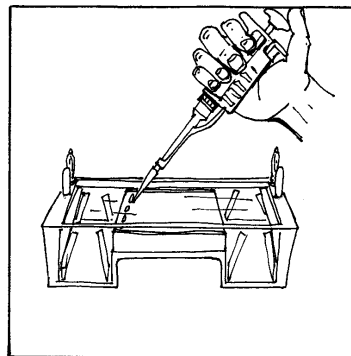
2. Connect the leads to the power supply, matching the color-coded red to red and black to black. **See Section 4.1 for recommended power conditions.**
3. Begin separation by electrophoresis.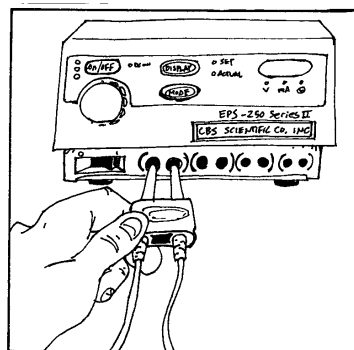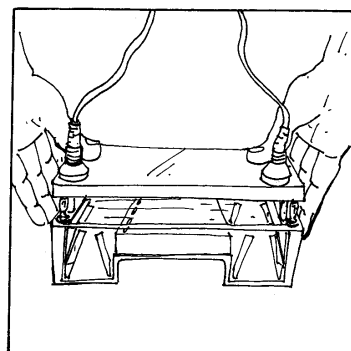

### 3.3 Removing the gel

Turn the power supply off and disconnect the leads from the power supply. Remove the safety cover from the Mini-Gel unit by placing thumbs on white posts next to red & black connectors, then pushing down while pulling up with fingers under lid. **DO NOT pull on power cords.**

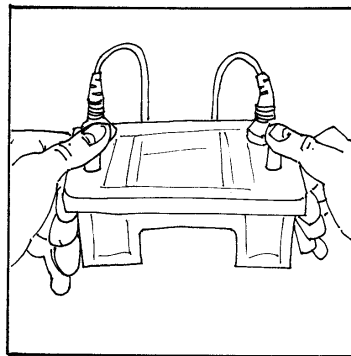Gently lift the gel or gel tray from the unit. **Always wear gloves, eye protection and protective clothing** if buffer and/or gel contain Ethidium Bromide. Ethidium Bromide is a powerful mutagen; gloves, eye protection and protective clothing should always be worn when handling the gel or buffer solutions. See Material Data Safety Sheets.

View separated fragments under UV light, using proper protection for eyes and skin (see manufacturer's instructions).

### 3.4 Maximum Well/Comb Volumes

NOTE: To calculate sample well volume expressed in millimeters (mm) of height, divide maximum volume by tooth depth.

MGC Comb for MGU-102T # of wells	Tooth width mm	Spacing between teeth mm	Overall length of comb mm	Tooth depth mm	1.0mm thickness volume per tooth microliters (ul)	2.0mm thickness volume per tooth microliters (ul)
1	49.4	1.57	49.4	6.35	313	616
3	15.4	1.57	49.4	6.35	97	195
8	4.77	1.57	49.4	6.35	30	60
10	3.5	1.57	49.4	6.35	22	44
12	2.6	1.57	49.4	6.35	16	32
14	2.0	1.57	49.4	6.35	12	24
16	1.6	1.57	49.4	6.35	10	20
MHC Comb for MGU-202T # of wells	Tooth width mm	Spacing between teeth mm	Overall length of comb mm	Tooth depth mm	1.0mm thickness volume per tooth microliters (ul)	2.0mm thickness volume per tooth microliters (ul)
1	69.8	1.5	69.8	6.35	443	886
3	22.2	1.5	69.8	6.35	141	282
5	12.6	1.5	69.8	6.35	80	160
8	7.3	1.5	69.8	6.35	46	92
10	5.5	1.5	69.8	6.35	34	68
12	4.3	1.5	69.8	6.35	27	54
14	3.5	1.5	69.8	6.35	22	44
16	2.8	1.5	69.8	6.35	17	34

## SECTION 4

### Running Conditions

#### 4.1 Recommended Power

The recommended power conditions for optimal resolution are 50 to 150V, constant, and 25 to 80 mA. The usual run time will vary for the voltage chosen but should range from 15 to 60 minutes, monitoring nucleic acid migration by progress of marker dyes. Constant power is not a necessity, but it produces uniform heat throughout the run. Be sure the polarity is correct i.e. that the DNA is loaded near the cathode (black electrode) to run toward the anode (red terminal).

Agarose gels may be stored for several days at 4°C wrapped in plastic wrap. Seakem Agarose (FMC) is used (normally) for preparative and analytical gels. Other types of agarose can be used for special purposes.

#### 4.2 Recommended Buffers

Type*	Concentrated Stock/liter	Final Concentration
TAE (Tris-acetate)	50X - 242 gm Tris base 57.1ml glacial acetic acid 100 ml 0.5M EDTA (pH 8.0)	1X - 0.04M Tris-acetate 0.001M EDTA
TBE (Tris-borate)	5X - 54 gm Tris base 27.5 gm boric acid 20 ml 0.5M EDTA (pH 8.0)	0.5X - 0.045M Tris-borate 0.001M EDTA

#### 10X - Loading Buffer (DNA)\*

0.25% Bromophenol blue  
0.25% Xylene cyanol  
20 % Ficoll Type 400  
0.1M EDTA, pH 8.0

### Ethidium Bromide Staining

One can add to the buffer only, the agarose gel solution only (after the solution has been made molten and the temperature has dropped to 55°C), or can be added after the electrophoresis run and the gel stained separately. Use at a final concentration of 0.1µg/ml from a 1mg/ml stock solution. Ethidium bromide is a Mutagen. ALWAYS wear gloves. Dispose of solutions in accordance with the safety regulations of your institution.

## 4.3 References

- Hames, B.D., Rickwood, D. (ed.) (1990). *Gel Electrophoresis of Nucleic Acids. A Practical Approach*. 2<sup>nd</sup> edn. IRL Press, Oxford. Ch 2.
- Sambrook, J., Fritsch, E.F., Maniatis, T. (1989). *Molecular Cloning. A Laboratory Manual*. 2<sup>nd</sup> edn. Cold Spring Harbor Laboratory Press, Cold Spring Harbor, New York. Ch 6.
- Ausubel, F.M., Brent, R., Kingston, R.E., Moore, D.D., Seidman, J.G., Smith, J.A., Struhl, K. (ed) (1993). *Current Protocols in Molecular Biology*. Vol. 1, Greene Publishing Associates, Inc. and John Wiley & Sons, Inc., Ch. 2.

## SECTION 5

### Maintenance of Equipment

#### 5.1 Care and Handling

The plastic components of the Mini-Horizontal units are fabricated from acrylic and polycarbonate. Electrodes and connectors are made from pure platinum, stainless steel, and chrome plated brass. As with any laboratory instrument, adequate care ensures consistent and reliable performance.

After each use, rinse buffer chamber, gel tray and combs with de-ionized water. Wipe dry with a soft cloth or paper towel, or allow to air dry. Whenever necessary, all components may be washed gently with water and a non-abrasive detergent, and rinsed and dried as above. *Never* use abrasive cleaners, glass cleaning sprays or scouring pads to clean the components, as these will damage the unit and components.

Additional precautions:

- Do not autoclave or dry-heat sterilize the apparatus or components.
- Do not expose the apparatus or components to phenol, acetone, benzene, halogenated hydrocarbon solvents or alcohols.
- Avoid prolonged exposure of the apparatus or components to UV light.
- Do NOT treat with diethylpyrocarbonate (DEPC)-treated water for extended periods at 37°C. A brief rinse with DEPC-water is sufficient after a thorough wash, followed by a quick rinse in 50% ethanol.

#### 5.2 Maintenance

The following inspection and maintenance procedures will help maintain the safety and reliable performance of the Mini-Horizontal systems. Replacement parts can be ordered by calling 1-858-755-4959 or by contacting your local distributor.

- Banana plugs and power cords should be inspected regularly. If the banana plugs become loose or do not feel friction tight replace the plugs or power cords.
- Should power cord assemblies (connectors, wire or shrouds) show any signs of wear or damage (e.g. cracks, nicks, abrasions, or melted insulation), replace them immediately.
- The platinum wire is secured to the banana jack by compression between a stainless washer and the jack nut. The nut/washer interface should be tight and free of corrosion.

## SECTION 6

### Horizontal Mini-Gel Kits and Accessories

Cat. #	Item
MGU-102T	<b>Horizontal Mini-Gel Kit-CE</b> , with tapered gates. Gel bed dimensions are 5.5cm x 8.5cm. Includes one comb, please specify comb cat. # from list below.
MGU-102T-FL	<b>Horizontal Mini-Gel Kit-CE with Flip Lid</b> , with tapered gates. Gel bed dimensions are 5.5cm x 8.5cm. Includes one comb, please specify comb cat. # from list below.
MGU-202T	<b>Horizontal Mini-Gel Kit-CE</b> , with tapered gates. Gel bed dimensions are 7.5cm x 10cm. Includes one comb, please specify comb cat. # from list below.
MGU-202T-FL	<b>Horizontal Mini-Gel Kit-CE with Flip Lid</b> , with tapered gates. Gel bed dimensions are 7.5cm x 10cm. Includes one comb, please specify comb cat. # from list below.

#### Color Coded Midi-Gel Option:

Available for the MGU-102T and MGU-202T models. To specify color desired, please add the following color code to the end of the catalog number: -Y (Yellow), -P (Purple), -B (Blue), -PK (Pink), or -G (Green). For example, to specify a green MGU-102T place order for catalog # MGU-102T-G.

### Mini-Gel Accessories

#### MGU-102T combs

(Available in Polycarbonate or Teflon-coated aluminum\*)

Cat. #	Comb dimensions
MGC-1001	1mm x 1 well
MGC-1003E	1mm x 3 equal wells
MGC-1003M	1mm x 2 well + middle marker
MGC-1008	1mm x 8 well
MGC-1010	1mm x 10 well
MGC-1012	1mm x 12 well
MGC-1014	1mm x 14 well
MGC-1016	1mm x 16 well
MGC-2001	2mm x 1 well
MGC-2003E	2mm x 3 equal wells
MGC-2003M	2mm x 2 well + middle marker
MGC-2008	2mm x 8 well
MGC-2010	2mm x 10 well
MGC-2012	2mm x 12 well
MGC-2014	2mm x 14 well
MGC-2016	2mm x 16 well

\* To specify Teflon combs add "-TF" to end of cat. #.

#### MGU-202T combs

(Available in Polycarbonate or Teflon-coated aluminum\*)

Cat. #	Comb dimensions
MHC-1001	1mm x 1 well
MHC-1003E	1mm x 3 equal wells
MHC-1003M	1mm x 2 well + middle marker
MHC-1008	1mm x 8 well
MHC-1010	1mm x 10 well
MHC-1012	1mm x 12 well
MHC-1014	1mm x 14 well
MHC-1016	1mm x 16 well
MHC-2001	2mm x 1 well
MHC-2003E	2mm x 3 equal wells
MHC-2003M	2mm x 2 well + middle marker
MHC-2008	2mm x 8 well
MHC-2010	2mm x 10 well
MHC-2012	2mm x 12 well
MHC-2014	2mm x 14 well
MHC-2016	2mm x 16 well

\* To specify Teflon combs add "-TF" to end of cat. #.

#### Dual Sided MGU-152T Combs (can be used with the MGU-102T)

(Available in Polycarbonate or Teflon-coated aluminum\*)

Cat. #	Comb dimensions
MGC-10-0810	1mm x 8 well/10 well
MGC-10-1216	1mm x 12 well/16 well
MGC-15-0810	1.5mm x 8 well/10 well
MGC-15-1216	1.5mm x 12 well/16 well
MGC-20-0810	2mm x 8 well/10 well
MGC-20-1216	2mm x 12 well/16 well

\* To specify Teflon combs add "-TF" to end of cat. #.

#### Dual Sided MGU-252T Combs (can be used with the MGU-202T)

(Available in Polycarbonate or Teflon-coated aluminum\*)

Cat. #	Comb dimensions
MHC-10-1014	1mm x 10 well/14 well
MHC-10-0816	1mm x 08 well/16 well
MHC-15-1014	1.5mm x 10 well/14 well
MHC-15-0816	1.5mm x 08 well/16 well
MHC-20-1014	2mm x 10 well/14 well
MHC-20-0816	2mm x 08 well/16 well

\* To specify Teflon combs add "-TF" to end of cat. #.

### Mini Gel Trays (UV Transparent)

Cat. #	Item
MGT-122	Gel Tray for the MGU-102T
MGT-222	Gel Tray for the MGU-202T

### Winged Gel Trays

Cat. #	Item
WGT-5585	Winged gel tray, 5.5cm x 8.5cm
WGT-7510	Winged gel tray, 7.5cm x 10cm

### Replacement Baffles

Cat. #	Item
BAFFLES-102T	Tapered Baffles for MGU-102T
BAFFLES-202T	Tapered Baffles for MGU-202T

### Gel Sealing Tape

Cat. #	Item
GT-72-10	1" thick x 72 yards
GT-72-15	1.5" thick x 72 yards

### Mini-Horizontal Gel Caster, MGU-102T

C.B.S. ☆ Scientific

### Mini-Horizontal Gel Caster, MGU-202T

Horizontal Mini-Gel

Cat. #	Item
GCT-101	Gel casting tray, single
GCT-103	Gel casting tray, three
GCT-106	Gel casting tray, six

Cat. #	Item
GCT-201	Gel casting tray, single
GCT-203	Gel casting tray, three
GCT-206	Gel casting tray, six

### Power Supply

Cat. #	Item
EPS-300-II	Mini Power Supply, with timer, CV or CC, 10-300V, 110V/60Hz, current range:4-500mA, 90 watts
EPS-300-IIV	Mini Power Supply, with timer, CV or CC, 10-300V, 220V/50Hz, current range:4-500mA, 90 watts
EPS-3000-III	Constant Power Supply, 0-3000 Volt, 110V/60Hz, Series II
EPS-3000-IIV	Constant Power Supply, 0-3000 Volt, 220V/50Hz, Series II

---

---

## NOTES

---

---

---

---

## NOTES

---

---

# CONTACT US

**Telephone:**  
Local or International  
858-755-4959  
Toll Free: 858-755-0733

**Fax:** 858-755-0733

**Check out our new and improved online ordering!**

[www.cbsscientific.com](http://www.cbsscientific.com)

**E-mail address:**  
[sales@cbssci.com](mailto:sales@cbssci.com)

**Mailing address:**  
C.B.S. Scientific Company  
P.O. Box 856  
Del Mar, CA 92014

**Shipping address:**  
C.B.S. Scientific Company  
10805 Vista Sorrento Pkwy  
Suite 100  
San Diego, CA 92121

**Credit Card Options:**  
Visa/Mastercard  
Discover/American Express